CONSORT

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

C731

October 1997

Index

Introduction	1
Manufacturer	1
Warranty	1
Servicing	1
Specifications	2
Keyboard	3
Error Codes	3
SET-UP	4
Channels	4
Date	4
Time	5
Temperature	5
Recorder	5
Alarm	6
Relay	6
Alternating measurements	7
Specific features	7
DATA ACQUISITION	7
Logging data	8
Adding data	8
Recalling data	8
Printing data	8
Diagnosis	9
Inputs	9
pH-MEASUREMENT	9
Maintenance	9
pH-measurement	10
mV-MEASUREMENT	11
Maintenance	11
Calibration	11
mV-measurement	11
CONDUCTIVITY MEASUREMENT	12
Maintenance	12
Calibration	12
Conductivity measurement	12
DATA COMMUNICATION	13
Recorder output	13
Digital output	13
Printer connection	13
Computer connection	14
TROUBLE SHOOTING	16
Simple test	16
Calibrate	16
Reset	16

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This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

Introduction

The normal operating temperature should be between 4° and 40°C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

Connect the instrument only to an earthed power line. The required power source is indicated on the label at the back of the instrument. Do not cut and splice the power cord. When removing the power cord from the wall outlet, be sure to unplug by holding the plug attachment and not by pulling the cord. Do not hold the plug by wet hand.

CONSORT nv
Parklaan 36
B2300 Turnhout
Belgium

Tel (++32)(14)41 12 79
Fax (++32)(14)42 91 79

Manufacturer

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

Warranty

CONSORT is not liable for consequential damages arising out of the use or handling of its products.

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply lead and **NOT** to send the following items unless they are suspect:

Servicing

Manual
Cables
Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

Specifications

Ranges	pH mV	0...14 pH ±2000 mV
	Conductivity	0...200 mS/cm
	Temperature	0...100°C
Resolution		0.01 pH, 1 mV, 0.1°C, 0.01 µS/cm
Inputs		2 BNC inputs for pH/mV/Ion electrodes 1 BNC input for a 2-pole conductivity cell 1 DIN input for a 4-pole conductivity cell 1 BANANA input for a Pt1000 temperature probe
Temp. Comp.		automatic with Pt1000 or manual
Digital output		programmable RS232, 150...4800 b/s, for bi-directional communication with a computer or printer
Analogue output		recorder-1: 0...2 V, ca 8 kΩ, programmable 8 bit D/A with adjustable span recorder-2: 0...2 V, ca 8 kΩ, programmable 8 bit D/A with adjustable span
Relays		1 programmable mechanical relay for high/low alarm, voltage-free contacts max. 250 V~/1A
Display		8+4 digit/character, 13 mm LED display, shows readings, °C & messages
Keys		8 tactile membrane keys
Ambient temp.		4...40°C
Rel. Humidity		0...90 % (non-condensing!)
Power supply		210-250 V~ (* 105-125 V~), 50/60 Hz, max. 4 VA
Cabinet		rugged plastic/metal cabinet
Dimensions		300 x 180 x 120 mm
Weight		2.4 kg

Keyboard

- MODE** = Selects all modes or escapes from error traps, calibration procedures, etc..by returning to the original mode.
- CAL** = Starts or proceeds a calibration or a function.
- ▲/▼** = Button for entering a value or for selecting a function.
- SET** = Installs the meter according your application.
- HOLD** = Holds display when measuring.
- RES** = Changes resolution/units of the display.
- PRINT** = Displayed value is printed through the RS232 output.
- ON/OFF** = Switches the instrument on or off.

Each time the instrument is switched on, it will check briefly its internal calibrations while showing all functions of the display and the software-version. An error message appears when something is wrong.

Error codes

- [Or. pH]** = Overrange pH (e.g. interrupted electrode cable).
- [Or. mV]** = Overrange mV (e.g. interrupted electrode cable).
- [Or. μ S]** = Overrange μ S (e.g. short circuited electrode cable).
- [Or. MS]** = Overrange mS (e.g. short circuited electrode cable).
- [Or. °C]** = Overrange °C (measured degrees out of range).
- [no dAtA]** = Data-logging memory empty.
- [not.CAL]** = Calibration procedure interrupted (the previous calibration remains valid!).
- [Err.SLP]** = Slope not within 80...120 % (e.g. worn pH electrode or used buffers).
- [Err.pHo]** = ISO pH not within 6...8 pH (e.g. defective electrode).
- [Err.cc]** = Cell constant not within 0.5...1.5 cm⁻¹.
- [Err.CAL]** = Calibration error (check electrodes and standard solutions).
- [Err.CH]** = Function not allowed for the selected channels.
- [Err.COd]** = Invalid code (enter the proper sequence of keys!).
- [Err.COM]** = General communication error.
- [Err.MEM]** = Fatal memory error (all factory settings are cleared. Have your instrument serviced!).
- [ILLEGAL]** = An unauthorised person has failed to repair the unit.
Warranty is elapsed from now on!

SET-UP

Channels

Procedure to pre-set the different channels and their possibilities according your applications.

[CH1=PH] : pH-measurement
[CH1=MV] : mV-measurement
[CH1=OFF] : Channel switched off
[EC=ON] / [EC=OFF] : Conductivity
[AtC=ON] / [AtC=OFF] : Automatic temperature compensation
[MtC=ON] / [MtC=OFF] : Manual temperature compensation

1. Press **SET** until the display shows [SEt CHAn] (choosing channels) and press **CAL** to continue (skipped when starting from a measuring mode).
2. The display shows [CH1=PH], [CH1=MV] or [CH1=OFF]. Select the desired mode for channel-1 with ▲ or ▼ and press **CAL** to continue or **SET** to return.
3. Go on in the same way with all next channels by repeating the procedure from point 2. Always press **CAL** to continue or **SET** to return.
4. The display shows [EC=on] or [EC=OFF]. Select the desired mode for the conductivity channel with ▲ or ▼ and press **CAL** to continue or **SET** to return.
5. This step is **skipped** unless the conductivity channel is on. The display shows [AtC=on], [MtC=on] or [AtC=OFF], [MtC=OFF]. Switch the temperature compensation on or off (for absolute measurements) with ▲ or ▼ and press **CAL** to continue or **SET** to return.
6. The display shows [rEF=25°C] or [rEF=20°C]. Select the temperature to which all future conductivity measurements will be referred with ▲ or ▼ and press **CAL** to continue or **SET** to return.

Procedure to change or display the date. When switched off, the date remains resident in the memory for logging and printing purposes.

Date

1. Press **SET** until the display shows [SEt dAtE] and press **CAL** to continue.
2. The display shows [dAtE=on] or [dAtE=OFF]. Select the desired mode with ▲ or ▼ and press **CAL** to continue or **SET** to return.
3. Program the date with ▲ or ▼ (respectively year, month and day) by pressing **CAL** each time (or press **SET** to return).

Time

Procedure to change or display the time. When switched off, the time remains resident in the memory for logging and printing purposes.

1. Press **SET** until the display shows [SEt tiME] and press **CAL** to continue.
2. The display shows [tiME=on] or [tiME=OFF]. Select the desired mode with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return.
3. Program the time with \blacktriangle or \blacktriangledown (respectively hour, minutes and seconds) by pressing **CAL** each time (or press **SET** to return).

This procedure permits to adjust the manual or automatic temperature compensation. For very accurate temperature readings the attached ATC-probe (Pt1000) can be calibrated in a $\pm 10^{\circ}\text{C}$ span.

Temperature

1. Press **SET** until the display shows [SEt $^{\circ}\text{C}$] and press **CAL** to continue.
2. When a Pt1000 temperature probe is connected to the instrument the display will show the measured temperature e.g. [$^{\circ}\text{C}$ =23.7][Aut]. If necessary, calibrate your probe with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return. Press **RES** to reset the temperature calibration to the original factory calibration.
3. When no Pt1000 temperature probe is connected to the instrument the display will show the manual temperature e.g. [$^{\circ}\text{C}$ =23.7][MAn]. Shift to the proper temperature with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return.

Two recorder outputs can be programmed individually to monitor the measurements of any channel.

Recorder

1. Press **SET** until the display shows [SEt rEC] and press **CAL** to continue.
2. The display shows e.g. [rc1=PH1], [rc1=MV2], [rc1=EC], [rc1= $^{\circ}\text{C}$], ...etc. Select the desired mode with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return.
3. This step is **skipped** unless [rc1=EC] has been chosen. The display shows the conductivity measuring range & units, e.g. [rc1 0.01MS][Unit]. Select the desired range with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return.
4. The display shows e.g. [rc1=6.00][LoW]. Select the desired minimum level corresponding to a 0 V recorder output with \blacktriangle or \blacktriangledown and press **CAL**.
5. The display shows e.g. [rc1=8.00][HiGH]. Select the desired maximum level corresponding to a 2 V recorder output with \blacktriangle or \blacktriangledown and press **CAL**.
6. The display shows e.g. [rc2=PH1], [rc2=MV2], [rc2=EC], [rc2= $^{\circ}\text{C}$], ...etc. Select the desired mode with \blacktriangle or \blacktriangledown and press **CAL** to continue in the same way for recorder-2 or **SET** to return. The display shows [SEt rEC].

Alarm

Procedure to pre-set high/low alarm levels for each channel. It will sound an alert or close a relay when readings stray outside the pre-set limits.

1. Press **SET** until the display shows [SEt ALrM] and press **CAL** to continue.
2. The display shows e.g. [ALrM PH1], [ALrM MV2], [ALrM EC], [ALrM °C]... etc. Select the desired alarm mode with ▲ or ▼ and press **CAL** to continue or **SET** to return.
3. The display shows [AL.=on] or [AL.=OFF]. Select the desired mode with ▲ or ▼ and press **CAL** to continue or **SET** to return.
4. This step is **skipped** unless [ALrM EC] has been chosen. The display shows the conductivity measuring range & units, e.g. [Lo 0.01MS][Unit]. Select the desired low-alarm range with ▲ or ▼ and press **CAL** to continue or **SET** to return.
5. The display shows e.g. [AL.=6.00][LoW]. Select the desired minimum level with ▲ or ▼ and press **CAL**. Each time the measurements are below this level, the beeper will sound an alarm (see 'Relay' page 6).
6. This step is **skipped** unless [ALrM EC] has been chosen. The display shows the conductivity measuring range & units, e.g. [Hi 0.01MS][Unit]. Select the desired high-alarm range with ▲ or ▼ and press **CAL** to continue or **SET** to return.
7. The display shows e.g. [AL.=8.00][HiGH]. Select the desired maximum level with ▲ or ▼ and press **CAL**. Each time the measurements are above this level, the beeper will sound an alarm (see 'Relay' page 6).

This instrument is provided with a mechanical relay. Its voltage-free contact can be programmed for high/low alarm or for instructing an automatic sampler to proceed when readings are stable.

Relay

1. Press **SET** until the display shows [SEt rEL] and press **CAL** to continue.
2. The display shows e.g. [rEL=ALrM] or [rEL=OFF]. Select the desired alarm mode with ▲ or ▼ and press **CAL** to continue or **SET** to return. By selecting [rEL=ALrM] the relay will close at each alarm.

Permits to install specific features such as beeper, baud-rate, printer-interval, identification-number and secret personal code.

1. Press **SET** until the display shows [SEtUP] and press **CAL** to continue.
2. The display lets you choose with ▲ or ▼ between [bEEP on] or [bEEP OFF]. Decide whether or not the beeper should sound each time the keyboard is touched. Press **CAL** to continue or press **SET** to return.
3. Press **CAL** until the display shows e.g. [br=2400]. Pre-set the desired baud-rate (150...4800 b/s) with ▲ or ▼ and press **CAL** to continue or press **SET** to return.
4. Press **CAL** until the display shows e.g. [rS=120]. Select the desired time interval (0...9999 s) between the transmitted data (RS232), with ▲ or ▼. Pre-set to zero if no automatic printing is required. Press **CAL** to continue or press **SET** to return.
5. Press **CAL** until the display shows e.g. [Id.no=5]. Select the desired identification number (0...999) for your meter with ▲ or ▼ and press **CAL** to continue or press **SET** to return. The instrument can be identified by e.g. a computer when a specific number is allocated to it. Pre-set to zero to ignore this number.
6. The display shows [Code on], [Code OFF] or [Code CAL]. Decide, with ▲ or ▼, whether all future access to the instrument should be by entering a code (on) or not (off). Press **CAL** to continue or press **SET** to return. Select [Code CAL] to enter your personal code to avoid undesired access to the instrument. The display shows [E=-----] while [=] is blinking. Enter your secret sequence of 5 keys to proceed. Press **CAL** to continue or **SET** to return.

DATA-ACQUISITION

Up to 500 values can be stored in the non-volatile memory of the instrument.

- [LOG dAtA]: Procedure to pre-set and start the data-logging.
[LSt dAtA] : Permits to view the stored values on the display.
[Prt dAtA] : Permits to send the stored values to a printer or computer.
[Cnt LOG] : Permits to continue with a next series of measurement.

[Log dAtA]: Procedure to pre-set and start the data-logging.

Logging data

1. Press **SET** until the display shows [SEt LoG] and press **CAL** to continue.
2. Select [Log dAtA] with \blacktriangle or \blacktriangledown and press **CAL** to continue or **SET** to return.
3. Choose with \blacktriangle or \blacktriangledown between [LOG PH1], [LOG PH2],..., [LOG μ S], ... etc. Select the desired mode and press **CAL**.
4. The display shows e.g. [dt=100]. Select the desired time interval (0...9999 s) between the data-logging, with \blacktriangle or \blacktriangledown and press **CAL**.
5. This step is **skipped** unless manual data-logging is chosen. Pre-set the time-interval (dt) to zero and press **CAL** to continue (point 6 is skipped).
6. The display shows e.g. [n=40]. Select the desired number of values (1...500) to be data-logged with \blacktriangle or \blacktriangledown and press **CAL**.
7. The display shows [StArt ?]. Press **CAL** to start the data-logging according the previous settings (a blinking [L] appears) or press **MODE** to return. After the last value has been data-logged the display will show [End LOG]. When manual data-logging has been selected, press **PRINT** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [LOG 7]. Press **SET** to return.

[Cnt LoG]: Procedure to continue with a next series of measurements, without erasing the previous data-logging.

Adding data

1. Press **SET** until the display shows [SEt LoG] and press **CAL** to continue.
2. Select [Cnt LoG] with \blacktriangle or \blacktriangledown provided the memory has enough capacity left! Press **CAL** to continue to continue with a next series of measurements, without erasing the previous data-logging.
3. Press **SET** to return.

[LSt dAtA]: Permits to view the stored values on the display.

Recalling data

1. Press **SET** until the display shows [SEt LoG] and press **CAL** to continue.
2. Select [LSt dAtA] with \blacktriangle or \blacktriangledown and press **CAL**.
3. The display shows the first data-logged value in memory (a blinking [n] appears). Press \blacktriangle or \blacktriangledown to view all other values.
4. While pressing **HOLD** the corresponding date is shown.
5. While pressing **RES** the corresponding time shown.
6. Press **PRINT** to send the recalled data to a printer.
7. Press **SET** to return.

[Prt dAtA]: Permits to send the stored values to a printer or computer.

Printing data

1. Press **SET** until the display shows [SEt LoG] and press **CAL** to continue.
2. Select [Prt dAtA] with \blacktriangle or \blacktriangledown and press **CAL** to print the data.
3. The display shows [PrintinG] while all data-logged values, including corresponding dates and times, are sent to the RS232 output in the following format, compatible with EXCEL™, LOTUS 123™,....The columns are separated by a TAB character:

Identification No: 037

#	VALUE	UNIT	C	CH	H	D
0001	7.34	pH	23.7	2	13:19:57	29/01/95
0002	7.21	pH	23.7	2	13:20:57	29/01/95
0003	7.08	pH	23.8	2	13:21:57	29/01/95
0004	6.71	pH	23.8	2	13:22:57	29/01/95
0005	6.49	pH	23.9	2	13:23:57	29/01/95

Permits to diagnose the performance of the electrodes.

Diagnosis

1. Press **SET** until the display shows [diAGn].
2. Sequentially press **CAL** to show the calibration details of all connected electrodes, e.g:

[94.6%] (slope)
[6.89 pHo] (ISO-potential)
[cc=1.034] (cell constant)

3. Press **CAL** to continue or press **SET** to return.

The pH/mV electrodes should be connected to the coaxial pH/mV connectors (1 & 2). If separate electrodes are used, connect the reference electrodes to the REF. terminals. Connect a 2-pole conductivity cell to the coaxial EC input or connect a 4-pole conductivity cell to the DIN input. **Never connect both types at the same time to the instrument!** Automatic temperature compensation and temperature measurements are possible by plugging a Pt1000 thermocompensator into the °C terminals. Without compensator, the manual temperature compensation is automatically switched on.

Inputs

pH-MEASUREMENT

A pH electrode is active and stable only after wetting! For this purpose it must be immersed for **at least ten hours** in a 3...4 M KCl solution. During short interruptions (e.g. storage) the electrode should be immersed in a 3...4 M KCl solution. In doing this it is always kept ready for use. When the interruption is longer than a month, refill the closing cap with 3...4 M KCl and plug it on the electrode tip in order to protect the glass bulb. Before use, ensure that the reference part of the electrode is topped up with a 3...4 M KCl solution.

Avoid a low pressure inside the electrode! Therefore always remove the closure from the refilling aperture during the measurements as well as during the standardisation. This allows the saltbridge solution to flow through the ceramic liquid junction and prevents contamination of the electrolyte. For the same reason, the inside level should always be higher than the outside level of the measuring solution. Close the refilling aperture again when storing the electrode.

A polluted electrode may be cleaned with a soft detergent or 0.1 M HCl. Greasy substances may be removed with acetone or alcohol (**never do this with plastic electrodes!**).

Maintenance

If the electrode is polluted by proteinaceous materials (such as blood), it should stand in a pepsin solution overnight and then be cleaned before use. The pH electrode wears away by being used. If the electrode tends to respond slower and calibration becomes difficult, even after cleaning, it should be replaced by a new one.

A pH measurement is always relative which makes it necessary to calibrate regularly.

pH measurement

Zero point:

Each electrode has its own specific pH value, corresponding to the zero potential point (ISO pH). This specific pH value can differ from its theoretical zero point (e.g. 7 pH) by a few tenths of a pH.

Slope:

A good electrode follows the Nernst equation within 90 to 100 %. If the slope lies outside these limits, the electrode becomes slow and may not be optimal for most applications.

1. Select the pH range by pressing **MODE** (see 'Channels' page 4). The display will immediately show the measured pH according to the previous standardisation. Should you want to restandardise, press **CAL**.
2. Rinse the electrodes with distilled water and immerse them in the first buffer solution.
3. The display shows one of the 9 buffers in memory (b1) or the manual input (M1) of user specified buffers, e.g. [b1=4.01] while [b1] or [M1] is blinking. Select the proper buffer with \blacktriangle or \blacktriangledown and press **CAL**. If you wish to change the manual buffer, first press **RES** to reset to the proper value with \blacktriangle or \blacktriangledown and then press **CAL**.
4. The instrument shows the measured buffer and will standardise automatically when readings are stable ([=] stops blinking). If the buffer has a different temperature, compensate the indicated value manually with \blacktriangle or \blacktriangledown (not necessary when using a Pt1000).
5. Go on in the same way with the next buffer (b2) and always press **CAL**. Press **MODE** to start the measurements.
6. After rinsing the electrodes with distilled water, immerse them in the samples and read the pH-value from the display.
7. Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.

- * See page 66 for the buffers in memory.
- * A blinking decimal point warns you for unstable measurements. Wait to read the display!
- * Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.
- * The instrument will refuse automatic standardisation when the electrode is unstable. Insufficient stirring or a worn electrode may be the cause. Select a lower resolution for a more rapid standardisation or replace the electrode!
- * Press **RES** to change the resolution from 0.1 to 0.01 pH.

mV-MEASUREMENT

Metal electrodes (Pt, Ag, Au): Metal electrodes are always ready for use. During short interruptions they are immersed in distilled water. **They should be cleaned regularly:**

- * Silver electrodes are immersed in a concentrated ammonia solution during one hour.
- * Platinum or gold electrodes are immersed in concentrated nitric acid during one hour.

1. Select the mV range by pressing **MODE** (see 'Channels' page 4).
2. After rinsing the electrodes with distilled water, immerse them in the solution to be measured and read the display.
3. Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.

Maintenance

***mV
Measurement***

CONDUCTIVITY MEASUREMENT

A conductivity cell is active and stable only after wetting! For this purpose it must be immersed for **at least one hour** in distilled water. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

A polluted cell may be cleaned with a soft detergent or diluted nitric acid. Greasy substances may be removed with acetone or alcohol (never do this with plastic electrodes!).

The cell constant depends on the physical dimensions of the platinum plates. They are coated with a spongy black platinum deposit (2-pole cell only!). This increases greatly the effective surface and reduces polarising effects. Since the coating of platinum black is quite resistant to chemical action, the cell constant remains fairly constant during the electrode life time.

1. Select the conductivity range by pressing **MODE** (the display shows [μ S] or [MS]) and press **CAL**.
2. The display will show the cell constant (e.g. [cc=1.073]). Press **CAL** to start the automatic determination procedure.
3. The display shows one of the KCl standard solutions in memory [St=0.01M] (1413 μ S/cm), [St=0.1M] (12.88 mS/cm), [St=1M] (111.8 mS/cm) or the manual input [St=MAN]. Select the proper standard with \blacktriangle or \blacktriangledown .
4. This step is **skipped** unless the manual input has been chosen. The display shows e.g. [MS=10.00] while the units are blinking. Choose with \blacktriangle or \blacktriangledown between 1.000 (units), 10.00 (tens) or 100.0 (hundreds) the desired calibration range and press **CAL** again. The display shows e.g. [MS=1.234] while [=] is blinking. Select the desired standard value with \blacktriangle or \blacktriangledown .
5. After rinsing the cell several times with the standard solution, immerse it the standard solution. The solution temperature is not so critical but it should lie between 0°C and 30°C. If the standard has a different temperature, compensate the indicated value manually with \blacktriangle or \blacktriangledown (see 'Temperature' page 5) (not necessary when using a Pt1000). Press **CAL** to continue.
6. The display will show the computed cell constant e.g. [cc=1.045] and will calibrate automatically when readings are stable ([=] stops blinking).

1. Select the conductivity range by pressing **MODE** (the display shows [μ S] or [MS]).
2. Rinse the electrodes with the sample and immerse them in it.
3. Read the display (remember: compensated to the reference temperature!). The autoranging feature gives you the largest possible number of digits + their unit.
4. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

Maintenance

Calibration

Conductivity measurement

- * See page 66 for the standards in memory.
- * A blinking decimal point warns you for unstable measurements. Wait to read the display!
- * Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.

DATA COMMUNICATION

Two recorders can be connected to both pairs red (+) and black (-) terminals. Use only laboratory recorders with a high input impedance! (See also 'Recorder' page 5).

Recorder output

A standard RS232 output terminal (DP9) is provided for interfacing the instrument with a printer or computer. Data is sent in the ASCII code at a BAUD rate of 150...4800 b/s (8 bit, no parity, 1 start and 2 stopbits, automatic line feed ON, data transfer protocol Xon/Xoff).

Digital output

Serial port pinout specifications:

- pin 1 : RLSD, received line signal detector
- pin 2 : TxD, transmit data
- pin 3 : RxD, receive data
- pin 4 : not connected
- pin 5 : Gnd, signal ground
- pin 6 : DSR, data set ready
- pin 7 : not connected
- pin 8 : CTS, clear to send
- pin 9 : not connected

Interfacing with a printer:

Press **PRINT** each time you want to print the measured values. The columns are separated by a TAB character. Run the procedure for automatic printing at timed intervals, 'Specific features' page 7, while you are measuring. The **PRINT** keypath does not function in this case. It is also disabled during an automatic data-logging., e.g.:

Printer connection

Identification No: 037

#	VALUE	UNIT	C	CH	H	D
0001	7.34	pH	23.7	2	13:19:57	29/01/95
0002	7.21	pH	23.7	2	13:20:57	29/01/95
0003	7.08	pH	23.8	2	13:21:57	29/01/95
0004	6.71	pH	23.8	2	13:22:57	29/01/95
0005	6.49	pH	23.9	2	13:23:57	29/01/95

Using WINDOWS™ to collect data:

1. Connect the computer with the RS232 terminal, using the optional AK2310 cable (eventually with adapter AB3190 for connection with a 25 pole RS232 terminal).
2. Start **WINDOWS™** and select *Terminal* in the *Accessories Group*.
3. Choose *Receive Text File* in the *Transfer Menu* and create a file name (TXT format) for the data to receive.
4. Select the *Settings Menu* and make sure the following parameters are programmed:

Baud Rate : 2400
Data bits : 8
Stop bits : 2
Parity : None
Flow control : Xon/Xoff

5. Start the instrument in the desired measuring mode for manual or automatic data collection (see 'Specific features' page 7 for more details).
6. When finished, open the created file with a standard spreadsheet (e.g. **EXCEL™**) or wordprocessor (e.g. **WORD™**) for further processing.

Interfacing with a computer:

Send a single character to the instrument and it will execute the corresponding command as follows:

- 1 = press **MODE**.
- 2 = press **CAL**.
- 3 = press **▲**.
- 4 = press **▼**.
- 5 = press **SET**.
- 6 = press **HOLD**.
- 7 = press **RES**.
- 8 = press **PRINT**.
- ? = send display to computer.
- + = keyboard on (manual operation remains possible).
- = keyboard off (manual operation disabled).
- V = start the input of a value (only valid in certain routines).

Follow exactly the same measuring or calibration procedures, as described in this manual, to include any desired command in your computer data-acquisition program.

This example continuously prints the measured values (pH and temperature) on the computer screen:

```
10 ON ERROR GOTO 70           'makes sure no error occurs
20 OPEN "COM1:2400,N,8,2" AS #1 'prepares computer input
30 PRINT #1,"8"              'PRINT button pressed
40 INPUT #1,V$               'measured values transmitted
50 PRINT V$                  'received values printed
60 GOTO 30                   'repeat procedure
70 RUN
```

Input of a value:

The direct input of values ("V" command) is only possible when the instrument is running a routine where also a manual input of these values could be made. If not, it will ignore any "V" command from the computer!

Examples of allowed routines are:

- * temperature input (e.g. manual temperature compensation).
- * control parameter input such as Low, High, ...
- * time settings.
- * input of manual buffer values.

Examples of NOT allowed routines are:

- * during measurements.
- * when values have to be selected rather than changed (e.g. buffer values in memory).
- * while an electrode is being calibrated.

A 16-bit value (2's complement) should be transmitted in the following sequence:

```
1st character = "V" (start a value input)
2nd character = highest byte of the value in ASCII
3rd character = lowest byte of the value in ASCII
4th character = checksum of 2nd and 3rd character in ASCII
5th character = LF, linefeed (ASCII-10)
```

When a correct checksum has been received, the instrument will send a confirmation character "!" to the computer. If not, a question mark "?" is sent. Both are eventually preceded by the identification number.

This example makes the computer to enter the value "1000" on instrument #7, in the sophisticated mode:

$$1000 = (\text{H-byte } 3) \times 256 + (\text{L-byte } 232)$$

```
10 OPEN "COM1:2400,N,8,2" AS #1 'prepares communication port
20 PRINT #1,"V";                'start a value input
30 PRINT #1,CHR$(3);            'highest byte of value
40 PRINT #1,CHR$(232);          'lowest byte of value
50 PRINT #1,CHR$(235);          'checksum of both bytes
60 PRINT #1,CHR$(10);           'LF, closes command
```

TROUBLE SHOOTING

Try this procedure if you suspect the electronic performance of the instrument.

Simple test

1. Short-circuit all pH/mV inputs. Leave all other inputs open.
2. Switch the instrument ON.
3. Select the conductivity range by pressing **MODE**. The display should show $[0.00\mu\text{S}] \pm 0.01 \mu\text{S/cm}$. Run the **Calibrate** procedure if erroneous values are read.
4. Select the mV ranges by pressing **MODE**. The display should show $[0\text{MV}] \pm 1 \text{mV}$. Run the **Calibrate** procedure if erroneous values are read.
5. Connect a fresh standard alkaline battery of 1.5 V to pH/mV input.
6. The display should now read approximately $[1500 \text{MV}]$. Run the **Reset** procedure if erroneous values are read.
7. Reverse the battery of 1.5 V and connect it to pH/mV input.
8. The display should now read approximately $[-1500 \text{MV}]$. Run the **Reset** procedure if erroneous values are read.

This procedure permits to calibrate automatically all amplifiers of the instrument.

Calibrate

1. Short-circuit all pH/mV inputs. Leave all other inputs open.
2. Switch the instrument ON while holding **CAL** pressed.
3. The display shows $[\text{SELFCAL}]$ for a while and returns to the measurements.

This procedure resets the instrument to the original factory settings. All personal memorised data (including access code) will be erased!

Reset

1. Switch the instrument ON while holding **MODE** pressed.
2. The display shows $[\text{rESEt}]$ for a while and returns to the measurements.

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