# CONSORT

---

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

C831

C832

C833

C835

C838

# QUALITYMASTERS

## CERTIFICATE OF REGISTRATION

This is to certify that:

### Consort NV

Parklaan 36  
B-2300 Turnhout

Has been assessed by QualityMasters in respect of its Quality Management System and found to comply with:

NEN-EN-ISO 9001 : 2000

Approval is hereby granted for registration providing the rules and conditions relating to certification are observed at all times.

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Lack of fulfillment of conditions as set forth in the certification regulations may render this certificate invalid. The use of the accreditation mark indicates accreditation in respect to the activities covered by the accreditation number EU9706010F.

# Q

# M

# REGISTRATION

# CERTIFICATE OF

Authorized signatory

*A. LEWIS*  
QualityMasters B.V.

This certificate remains the property of QualityMasters B.V.

# WARRANTY CERTIFICATE

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory.

**CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse.

A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

## DECLARATION OF CONFORMITY

We declare under our sole responsibility that the product

**Multi-channel analysers**  
**content of the type numbers**  
**C831, C832, C833, C835, C838**

to which this declaration relates is in conformity  
with the following standards

**EN61010**  
**LOW VOLTAGE DIRECTIVE 73/23/EEG**  
**EN50081-1**  
**EN50082-1**  
**EN60555-2**  
**EMC DIRECTIVE 89/336/EEG**

Turnhout, November 10, 1995

A handwritten signature in black ink, appearing to be 'J. G. G.' or similar, is written over a horizontal line.

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**Forms**

Buffers, standards .....	69
Technical assistance form - pH .....	70
Technical assistance form - Conductivity .....	71
Technical assistance form - Ion.....	72
Technical assistance form - Oxygen .....	73
Your satisfaction .....	74

This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 4° and 40° C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

Connect the jack of a suitable AC adapter (model **A4070** for 230 V~ or model **A4071** for 115 V~) to the DC socket. Do not hold the adapter by wet hand.

---

**Manufacturer****CONSORT nv**

Parklaan 36  
B2300 Turnhout  
Belgium

Tel (+32)(14)41 12 79  
Fax (+32)(14)42 91 79  
E-mail: info@consort.be

---

**Warranty**

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

**CONSORT** is not liable for consequential damages arising out of the use or handling of its products.

---

**Servicing**

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply lead and **NOT** to send the following items unless they are suspect:

*Manual*

*Cables*

*Accessories*

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

<b>Ranges</b>	<p>pH -2...+16 pH</p> <p>mV <math>\pm 2000</math> mV</p> <p>Ion <math>10^{-11}</math> to <math>10^{-1}</math>, any unit (C833, C835)</p> <p>rH<sub>2</sub> 0...42 rH<sub>2</sub> (C838)</p> <p><math>\mu</math>W 0...400000 <math>\mu</math>W (C838)</p> <p><b>Conductivity</b></p> <p>cc = 0.1 cm<sup>-1</sup> : 0.001 <math>\mu</math>S/cm ... 20 mS/cm (not C831)</p> <p>cc = 1 cm<sup>-1</sup> : 0.01 <math>\mu</math>S/cm ... 200 mS/cm</p> <p>cc = 10 cm<sup>-1</sup> : 0.1 <math>\mu</math>S/cm ... 2000 mS/cm (not C831)</p> <p><b>Resistivity (C832, C833, C835, C838)</b></p> <p>cc = 0.1 cm<sup>-1</sup> : 100 <math>\Omega</math>.cm ... 200 M<math>\Omega</math>.cm</p> <p>cc = 1 cm<sup>-1</sup> : 10 <math>\Omega</math>.cm ... 20 M<math>\Omega</math>.cm</p> <p>cc = 10 cm<sup>-1</sup> : 1 <math>\Omega</math>.cm ... 2 M<math>\Omega</math>.cm</p> <p><b>TDS (C832, C833, C835, C838)</b></p> <p>cc = 0.1 cm<sup>-1</sup> : 0.01 mg/l ... 20 g/l</p> <p>cc = 1 cm<sup>-1</sup> : 0.01 mg/l ... 100 g/l</p> <p>cc = 10 cm<sup>-1</sup> : 0.1 mg/l ... 100 g/l</p> <p><b>Salinity</b> 0 ... 70 ppt (C832, C833, C835, C838)</p> <p><math>^{\circ}</math>C 0 ... 100<math>^{\circ}</math>C (O<sub>2</sub>: 0...50<math>^{\circ}</math>C)</p> <p>O<sub>2</sub> 0 ... 60 mg/l (C835)</p> <p><b>Saturation</b> 0 ... 600 % (C835)</p>
<b>Resolution</b>	0.01/0.001 pH, 0.1 mV, 0.01 rH <sub>2</sub> , 0.1 $^{\circ}$ C, 0.01/0.001 $\mu$ S/cm, 1/10 $\Omega$ .cm, 1% ion, 0.01 mg/l, 0.1 % O <sub>2</sub>
<b>Inputs</b>	<p>C831-C833: 2 BNC inputs for pH/mV/Ion electrodes</p> <p>C835: 6 BNC inputs for pH/mV/Ion electrodes (channel-2 = input for oxygen electrode)</p> <p>C838: 1 BNC input for pH, 1 BNC input for mV 1 BNC input for a conductivity cell 1 BANANA input for a Pt1000 temperature probe</p>
<b>Temp. Comp.</b>	automatic with Pt1000 or manual
<b>Digital output</b>	programmable RS232, 150...4800 b/s, for bi-directional communication with a computer or printer
<b>Analogue output</b>	0...2 V, ca 8 k $\Omega$ , programmable 8 bit DAC
<b>Display</b>	8+4 digit/character, 13 mm LED display
<b>Keys</b>	9 tactile membrane keys
<b>Ambient temp.</b>	4...40 $^{\circ}$ C
<b>Rel. Humidity</b>	0...90 % (non-condensing!)
<b>Power supply</b>	210-250 V- (* 105-125 V-), 50/60 Hz, max. 4 VA
<b>Cabinet</b>	IP65 cabinet
<b>Dimensions</b>	260 x 180 x 90 mm
<b>Weight</b>	1 kg

---

MODE	=	Selects all modes or escapes from error traps, calibration procedures, etc..by returning to the original mode.
CAL	=	Starts or proceeds a calibration or a function.
▲/▼	=	Button for entering a value or for selecting a function.
SET	=	Installs the meter according your application.
HOLD	=	Holds display when measuring.
RES	=	Changes resolution/units of the display.
PRINT	=	Displayed value is printed through the RS232 output.
ON/OFF	=	Switches the instrument on or off.

---

**Error codes**

[Or. pH]	=	overrange pH
[Or. mV]	=	overrange mV
[Or. µS]	=	overrange µS
[Or. MS]	=	overrange mS
[Or. OhM]	=	overrange Ω
[Or. Cn]	=	overrange concentration
[Or. °C]	=	overrange °C
[Or.PPt]	=	overrange ppt (TDS)
[Or.SAL]	=	overrange salinity
Or.%]	=	overrange %
[no dAtA]	=	data-logging memory empty
[not.CAL]	=	calibration procedure interrupted
[Err.SLP]	=	slope not within 80...120 %
[Err.pHo]	=	ISO pH not within 6...8 pH
[Err.cc]	=	cell constant not within 70...130% of the nominal value
[Err.CAL]	=	calibration error
[Err.Stb]	=	electrode not stable
[not drY]	=	cell too wet for automatic capacitive compensation
[diLutE]	=	solution too concentrated for a TC plotting
[Err.EC]	=	no cell connected during a TC plotting
[Err. °C]	=	initial temperature too high (>10°C) or no temperature probe connected
[Err.SAL]	=	salinity measurement error during O <sub>2</sub> -measurements
[Err.tC]	=	unprogrammed temperature-curve
[Err.Cn]	=	unprogrammed concentration-curve
[Err.COd]	=	invalid code
[Err.COM]	=	general communication error
[Err.MEM]	=	fatal memory error; have your instrument serviced!
[ILLEGAL]	=	an unauthorised person has failed to repair the unit; <b>warranty is elapsed from now on!</b>

---

**Inputs**

The pH/mV/Ion electrodes should be connected to the coaxial pH/mV connectors CH.1, etc. If separate electrodes are used, connect the reference electrodes to the REF. terminals. Connect the conductivity cell to the coaxial EC input. Automatic temperature compensation and temperature measurements are possible by plugging a Pt1000 thermocompensator into the °C terminals. Without compensator, the manual temperature compensation is automatically switched on. C835: the O<sub>2</sub>-electrode should be connected to the coaxial connector CH.2. Its banana plugs should be inserted in the °C terminals.

---

**Recorder output**

A recorder can be connected to the red (+) and black (-) terminals. Use only laboratory recorders with a high input impedance!

A standard RS232 output terminal (DP9) is provided for interfacing the instrument with a printer or computer. Data is sent in the ASCII code at a BAUD rate of 150...4800 bps (8 bit, no parity, 1 stopbit).

Serial port pinout specifications:

- pin 1 : connected to pin 4 and pin 6
- pin 2 : TxD, transmit data
- pin 3 : RxD, receive data
- pin 4 : connected to pin 1 and pin 6
- pin 5 : Gnd, signal ground
- pin 6 : connected to pin 1 and pin 4
- pin 7 : connected to pin 8
- pin 8 : connected to pin 7
- pin 9 : not connected

---

*(C831, C832, C833, C835)*

*Channels*

1. Press **SET** until the display shows [SEt CHAn] and press **CAL**.
2. The display shows [CH1=PH], [CH1=MV], [CH1=Cn] or [CH1=OFF]. Select the desired mode for channel-1 and press **CAL**.
3. The display shows [CH2=PH], [CH2=MV], [CH2=Cn], [CH2=MG.L] [CH2=%] or [CH2=OFF]. Select the desired mode for channel-2 and press **CAL**.
4. Go on in the same way with all next channels by repeating the procedure from point 2. Always press **CAL**.
5. **C832, C833**: this step is **skipped** unless channel-1 and channel-2 have been programmed for the same mode (pH or mV). The display shows [dpH=on] or [dmV=on], [dpH=OFF] or [dmV=OFF]. Select the desired mode and press **CAL**. Mode dpH/dmV permits to read the difference between both channels.
6. The display shows [EC=on] or [EC=OFF]. Select the desired mode for the conductivity channel and press **CAL**.
7. This step is **skipped** unless the conductivity channel is on. The display shows [AtC=on], [MtC=on] or [AtC=OFF], [MtC=OFF]. Switch the temperature compensation on or off (for absolute measurements) and press **CAL**.
8. The display shows [rEF=25°C] or [rEF=20°C]. Select the temperature to which all future conductivity measurements will be referred and press **CAL**.

---

*(C838)*

*Reference*

1. Press **SET** until the display shows [SEt rEF] and press **CAL** to enter the reference type of the redox electrode [CALOMEL] or [AgCL]. Press **CAL** to continue or **SET** to return.
2. The display shows e.g. [rEF=25°C]. Select the temperature (20...40°C) to which all future conductivity measurements will be referred and press **CAL**.

---

*Date*

1. Press **SET** until the display shows [SEt dAtE] and press **CAL**.
2. The display shows [dAtE=on] or [dAtE=OFF]. Select the desired mode and press **CAL**.
3. Program the date (respectively year, month and day) by pressing **CAL** each time.

---

*Time*

1. Press **SET** until the display shows [SEt tiME] and press **CAL**.
2. The display shows [tiME=on] or [tiME=OFF]. Select the desired mode and press **CAL**.
3. Program the time (respectively hour, minutes and seconds) by pressing **CAL** each time.

1. Press **SET** until the display shows [SEt °C]. Select the desired mode and press **CAL**.
2. The display will show the measured temperature e.g. [°C=23.7][Aut]. If necessary, calibrate your probe and press **CAL**. Press **RES** to reset the temperature calibration to the original factory calibration.
3. When no Pt1000 temperature probe is connected to the instrument the display will show the manual temperature e.g. [°C=23.7][MAn]. Shift to the proper temperature and press **CAL**.

## Recorder

1. Press **SET** until the display shows [SEt rEc] and press **CAL**.
2. The display shows for all channels e.g. [rEc=PH1], [rEc=MV2], [rEc=Cn1], [rEc=Cn2], [rEc=EC], [rEc=dPH], [rEc=MG.L], [rEc=%], [rEc=°C]...etc. Select the desired mode and press **CAL**.
3. This step is **skipped** unless [rEc=EC] has been chosen. Press **CAL** and the display shows the conductivity measuring range and units, e.g. [rEc 0.01mS][Unit]. Select the desired range and press **CAL**.
4. The display shows e.g. [rEc=6.00][LoW]. Select the desired minimum level corresponding to a 0 V recorder output and press **CAL**.
5. The display shows e.g. [rEc=8.00][HiGH]. Select the desired maximum level corresponding to a 2 V recorder output and press **CAL**. The display shows [SEt rEc].

## Specific features

1. Press **SET** until the display shows [SEtUP] and press **CAL**.
2. The display lets you choose between [bEEP on] or [bEEP OFF]. Decide whether or not the beeper should sound each time the keyboard is touched and press **CAL**.
3. The display lets you choose between [POWER][on] or [POWER][OFF]. Choose [POWER][on] to inactivate the ON/OFF-key (meter is always on and cannot be switched off). After an interruption of the mains power, the meter will resume measurements automatically. Choose [POWER][OFF] to activate the ON/OFF-key (meter can be switched on or off). Press **CAL**.
4. The display shows e.g. [br=2400]. Pre-set the desired baud-rate (300...9600 b/s) and press **CAL**.
5. The display shows e.g. [rS=120]. Select the desired time interval (0...9999 s) between the transmitted data (RS232). Pre-set to zero if no automatic printing is required and press **CAL**.
6. The display shows e.g. [Id.no=5]. Select the desired identification number (0...999) for your meter and press **CAL**. The instrument can be identified by e.g. a computer when a specific number is allocated to it. Pre-set to zero to ignore this number.
7. The display shows [Code on], [Code OFF] or [Code CAL]. Decide whether all future access to the instrument should be by entering a code (on) or not (off). Press **CAL**. Select [Code CAL] to enter your personal code to avoid undesired access to the instrument. The display shows [E=-----] while [=] is blinking. Enter your secret sequence of 5 keys to proceed. Press **CAL**.

## Alternating measurements

1. Press **SET** until the display shows [ALt] and press **MODE** to start the alternating mode. The display alternates between the selected modes with an interval of 4 s. Pressing **PRINT** in this mode will print all values of the selected channels.
2. Press **MODE** again to stop this mode and continue with the actual channel.

1. Select the pH range by pressing **MODE**. The display will immediately show the measured pH according to the previous standardisation. Should you want to recalibrate, press **CAL**.
2. **C833, C835**: choose between [Cor OFF], [Cor on] or [Cor CAL]. Decide whether or not temperature corrected readings should be displayed (normally OFF), press **CAL** and proceed with point 4. Select [Cor CAL] to plot the temperature curve of your sample in the memory of the instrument and press **CAL**.
3. **C833, C835**: the display shows e.g. [rEF=25°C] while [=] is blinking. Select the temperature to which all future measurements will be referred to and press **CAL** to start the automatic plotting procedure or press **MODE** to cancel. Prepare the sample solution to be measured in a thermostatic bath with magnetic stirrer and immerse the electrodes in it. Its temperature should be about 5°C lower than the lowest value you may possibly need during future measurements. The display shows e.g. [7.87cPH][18.9] which means that the sample has 7.87 pH at 18.9°C. A blinking [c], indicating that plotting has started, appears. Wait a while to enable the electrodes to adapt themselves to the new temperature. Press **HOLD**, when readings are stable, to store the values in memory. Increase the temperature with at least 5°C and repeat the same procedure for as many temperatures as necessary. When the temperature has reached about 5°C more than the highest value you need for future measurements, press **CAL** to stop the temperature tracing. The display shows briefly [WAlt] while the curve is being computed.
4. Rinse the electrodes with distilled water and immerse them the first buffer solution.
5. The display shows one of the 9 buffers in memory (b1) or the manual input (M1) of user specified buffers, e.g. [b1=4.01] while [b1] or [M1] is blinking. Select the proper buffer and press **CAL**. If you wish to change the manual buffer, first press **RES** to reset to the proper value and then press **CAL**.
6. The instrument shows the measured buffer and will standardise automatically when readings are stable ([=] stops blinking). If the buffer has a different temperature, compensate the indicated value manually (not necessary when using a Pt1000).
7. Go on in the same way with all the next buffers (b2...b5) and always press **CAL** or press **MODE** to start the measurements immediately.
8. After rinsing the electrode with distilled water, immerse them in the samples and read the pH-value from the display.
9. Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.

#### Relative pH measurements (not C838):

1. While in the pH range, press  $\blacktriangle$ . At that moment the instrument stores the actual pH and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
  2. To cancel relative pH readings, press **MODE** and you will be in the normal measuring mode again.
- *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
  - *Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.*
  - *Press SET to change the resolution from 0.1 to 0.001 pH.*

**Maintenance:**

A pH electrode is active and stable only after wetting! For this purpose it must be immersed for at least ten hours in a 3...4 M KCl solution. During short interruptions (e.g. storage) the electrode should be immersed in a 3...4 M KCl solution. In doing this it is always kept ready for use. When the interruption is longer than a month, refill the closing cap with 3...4 M KCl and plug it on the electrode tip in order to protect the glass bulb. Before use, ensure that the reference part of the electrode is topped up with a 3...4 M KCl solution.

Avoid a low pressure inside a refillable electrode! Therefore always remove the closure from the refilling aperture during the measurements as well as during the calibration. This allows the saltbridge solution to flow through the ceramic liquid junction and prevents contamination of the electrolyte. For the same reason, the inside level should always be higher than the outside level of the measuring solution. Close the refilling aperture again when storing the electrode.

A polluted electrode may be cleaned with a soft detergent or 0.1 M HCl. Greasy substances may be removed with acetone or alcohol (**never do this with plastic electrodes!**).

If the electrode is polluted by proteinaceous materials (such as blood), it should stand in a cleaning solution overnight and then be cleaned with distilled water before use. The pH electrode wears away by being used. If the electrode tends to respond slower and calibration becomes difficult, even after cleaning, it should be replaced by a new one.

---

**mV measurement**

1. Select the mV range by pressing **MODE**.
2. After rinsing the electrodes with distilled water, immerse them in the solution to be measured and read the display.
3. Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.

**Calibration:**

1. Select the mV range by pressing **MODE**.
2. Press **CAL** to adjust the mV scale with any reference. Immerse the electrodes in a standard solution of known potential.
3. The display shows e.g. [MV=476.3]. Adjust to the proper value and press **CAL** again or press **RES** to reset the absolute readings.

**Relative mV measurements (not C838):**

1. While in the mV range, press  $\blacktriangle$ . At that moment the instrument stores the actual mV and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
2. To cancel relative mV readings, press **MODE** and you will be in the normal measuring mode again.

- Press **RES** to change the resolution from 1 to 0.1 mV. (not C831)
- C838: press  $\blacktriangle$  or  $\blacktriangledown$  to switch between mV readings referred to a standard hydrogen electrode [MV.H] or not [MV].

**(C838)**

---

**rH<sub>2</sub> measurement**

1. Select the rH<sub>2</sub> range by pressing **MODE**.
2. After calibrating and rinsing the pH and redox electrodes with distilled water, immerse them in the solution to be measured and read the display.

1. Select the conductivity range by pressing **MODE** (the display shows [ $\mu$ S] or [MS]).
2. After rinsing the electrode with distilled water and then with some of the sample, immerse it the sample.
3. Read the display (remember: compensated to the reference temperature!). The autoranging feature gives you the largest possible number of digits + their unit.
4. Rinse the electrode always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

#### Calibration:

1. Select the conductivity range by pressing **MODE** (the display shows [ $\mu$ S] or [MS]) and press **CAL**.
2. The display will show the cell constant (e.g. [cc=1.073]). Press **CAL** to start the automatic determination procedure.
3. The display shows one of the KCl standard solutions in memory [St=0.01M][cc1] (1413  $\mu$ S/cm), [St=0.1M][cc1] (12.88 mS/cm), [St=1M][cc1] (111.8 mS/cm) or the manual input [St=MAN]. Select the proper standard.
4. This step is **skipped** unless the manual input has been chosen. The display shows e.g. [MS=10.00][cc1] while the units are blinking. Choose between 1.000 (units), 10.00 (tens) or 100.0 (hundreds) the desired calibration range and press **CAL** again. The display shows e.g. [MS=1.234][cc1] while [=] is blinking. Select the desired standard value (at real temperature!) (cc2 and cc3 should be at least 5 times stronger than the previous one).
5. After rinsing the electrode with distilled water and then with some of the standard solution, immerse it the standard solution. The solution temperature is not so critical but it should lie between 0°C and 30°C. If the standard has a different temperature, compensate the indicated value manually (not necessary when using a Pt1000). Press **CAL**.
6. The display will show the computed cell constant e.g. [cc1=1.045] and will calibrate automatically when readings are stable ([=] stops blinking).
7. Go on in the same way from step 5 with all the next standard solutions (cc2...cc3) by pressing **CAL** each time. Press **MODE** to stop the calibration procedure.

#### Capacitive compensation:

1. Capacitive compensation increases accuracy in the very low conductivity ranges (<10  $\mu$ S/cm). Verify if the attached electrode is completely dry.
  2. Select the conductivity range by pressing **MODE** (the display shows [ $\mu$ S] or [MS]) and press **CAL**.
  3. The display will show the cell constant (e.g. [cc=1.073]), [COMP.on] or [COMP.OFF]. Select [COMP.on] and press **CAL** to start the automatic capacitive compensation. The display will show [WaiT] for a while.
- *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
  - *C832, C833, C835, C838: press RES to switch between ppm (TDS: ppm = mg/l, ppt = g/l), SAL (salinity), S/cm (conductivity) or  $\Omega$ .cm (resistivity: E6 = M $\Omega$ , E3 = k $\Omega$ , E0 =  $\Omega$ ).*
  - *C832, C833, C835: when starting a titration, press  $\nabla$  to lock the actual range and avoid cross-over errors due to differences in the measuring frequency. Press MODE to return to the normal measuring mode.*
  - *Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.*

### Relative measurements (not C838):

1. While in the conductivity or salinity mode, press  $\Delta$ . At that moment the instrument stores the actual value and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
2. To cancel relative readings, press **MODE** and you will be in the normal measuring mode again.

### Maintenance:

A conductivity cell is active and stable only after wetting! For this purpose it must be immersed for at least one hour in distilled water. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

A polluted cell may be cleaned with a soft detergent or diluted nitric acid. Greasy substances may be removed with acetone or alcohol (never do this with plastic electrodes!).

### Temperature compensation (C833, C835):

*Each solution has its own temperature coefficient, which also varies with the temperature. Therefore it is impossible to achieve an exact temperature compensation over a wide span of temperatures with a conventional conductometer. However, this model features a unique microprocessor controlled system, which helps you to completely overcome the problem of non linear temperature compensation curves for most of the species to be measured. There is no need to keep sample temperatures close to each other, since you can plot special temperature curves from up to 10 different solutions.*

1. Press **MODE** until the display will show e.g. [tc Std] or [tc MAn]. Remember that [tc Std] is the fixed standard temperature curve (EN27888) which can never be altered. Select the desired temperature curve and press **MODE** to proceed or **CAL** to start the automatic tracing.
  2. Prepare the sample solution to be measured in a thermostatic bath with magnetic stirrer. Its temperature should be about 5°C lower than the lowest value you may possibly need during future measurements, but at least lower than 10°C. The display shows e.g. [tc MAn][8.9] which means manual temperature curve at 8.9°C. After rinsing the electrodes with the solution to be measured, immerse them in that solution. Wait a while to enable the electrodes to adapt themselves to the low temperature. Press **CAL** again to start the automatic plotting procedure.
  3. Slowly increase the temperature (max. 0.1°C/s) of the thermostatic bath. You can monitor the temperature on the display while a blinking [c], indicating that plotting has started, appears. When the temperature is about 5°C higher than the highest value you need for future measurements (min. 40°C required), press **CAL** to stop the temperature tracing. The display shows shortly [WAlt] while the curve is being computed.
- After a temperature curve has been selected and **PRINT** is pressed, the instrument transmits the whole curve to the digital output. In the meantime the display shows shortly [PrintinG].
  - If you have traced the temperature compensation in a limited range (e.g. from 10° to 40°C), the curve is supposed to go on linearly outside these limits.

(C838)

$\mu$ W measurement

1. Select the  $\mu$ W range by pressing **MODE**.
2. After calibrating and rinsing the pH, ORP and conductivity electrodes with distilled water, immerse them in the solution to be measured and read the power in  $\mu$ W.

**Calibration:**

Always calibrate the electrode in the same conditions as future determinations will require and make sure temperature remains constant. The instrument only accepts the following standard values, in scientific notation:  $1 \cdot 10^9$  /  $3 \cdot 10^9$  /  $1 \cdot 10^8$  /  $3 \cdot 10^8$  /  $1 \cdot 10^7$  /  $3 \cdot 10^7$  /  $1 \cdot 10^6$  /  $3 \cdot 10^6$  /  $1 \cdot 10^5$  /  $3 \cdot 10^5$  /  $1 \cdot 10^4$  /  $3 \cdot 10^4$  /  $1 \cdot 10^3$  /  $3 \cdot 10^3$  /  $1 \cdot 10^2$  /  $3 \cdot 10^2$  /  $1 \cdot 10^1$  /  $3 \cdot 10^1$  /  $1 \cdot 10^0$ , any unit. Therefore, e.g. 1 mg/l is shown on the display as [1.-3] (=  $1 \cdot 10^3$ ).

1. Prepare 2 or more (up to 19) standards, covering your future measuring range. Each of them should be 3 to 10 times stronger than the previous one.
  2. Press **MODE** until the display shows [Dir.Conc1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Dir.Conc1]. Press **CAL** to start the automatic plotting procedure.
  3. After rinsing the electrodes with distilled water, immerse them in the first standard solution. The display will show e.g. [3.-7 263.8] (=  $3 \cdot 10^7$  / 263.8 mV). Select the first calibration value and press **CAL** when readings are stable.
  4. After rinsing the electrodes with distilled water, immerse them in the second standard solution. The display will show e.g. [1.-6 206.5] (=  $1 \cdot 10^6$  / 206.5 mV). Select your next standard and press **CAL** when readings are stable.
  5. Go on in the same way with all the next standard solutions. After the last one has been entered, press **MODE** to stop the calibration procedure. The instrument will now compute all values to a curve. In the mean time the display shows briefly [WAlt].
  6. After rinsing the electrodes with distilled water, immerse them in a blank solution. Select [bL=on] or [bL=OFF] and press **CAL**. The display will show e.g. [Eo=85.3]. When readings are stable press **CAL**.
  7. The display shows one of the following modes: [Dir.Conc1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1]. Recall the desired method and press **MODE** to proceed.
- After a method has been selected and **PRINT** is pressed, the instrument transmits the whole curve to the digital output. In the meantime the display shows shortly [PrintinG].
  - Pressing **RES** during the measuring phase, permits you to choose between readings in scientific notation or in ng/l, µg/l, mg/l or g/l.

**Direct concentration:**

1. Press **MODE** until the display shows [Dir.Conc1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Dir.Conc1]. Press **MODE** to proceed.
2. Rinse the electrodes with distilled water, immerse them in the samples, and read the concentration on the display.

**Standard addition:**

*The standard addition method is used when the samples can be kept at the same temperature and a standard of 5 to 100 times the expected sample concentration is available. The most suitable increment can be found from: increment percentage > 100 x expected sample value / standard value.*

1. Press **MODE** until the display shows [Dir.CONC1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Std.Add1]. Press **MODE** to proceed.
2. The display shows one of the possible increments (1, 2, 5, 10, or 20 %), e.g. [10% Incr]. Select the desired value and press **CAL**.
3. Prepare a standard solution of the measured species. Its value should be 5 to 100 times higher than the expected sample concentration. The display shows one of the possible standards e.g. [Std 1.-3]. Select the desired value and press **CAL**.
4. After rinsing the electrodes with distilled water, immerse them in e.g. 100 ml sample solution. The display will show the electrode potential e.g. [E=307.6]. When readings are stable press **CAL**.
5. Add a small increment corresponding to the chosen mode (e.g. 10 ml). The display will show the difference in electrode potential e.g. [dE=48.1]. When readings are stable press **CAL**. The instrument calculates the result and shows the concentration of the sample in the same unit as the standard used for the addition.
6. For a next measurement press **CAL** and the same procedure starts again from point 4). Press **MODE** to abort.

**Standard subtraction:**

*The standard subtraction method is used to measure species for which stable standards or ion selective electrodes do not exist. The procedure is similar to the standard addition method. The most suitable increment can be found from: increment percentage > 100 x expected sample value / standard value.*

1. Calculate the stoichiometric ratio (n) of the reaction:  $xS + yR \rightarrow S_xR_y$   $n=x/y$  (S = sample, R = reagent).
2. Press **MODE** until the display shows [Dir.CONC1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Std.Sub1]. Press **MODE** to proceed.
3. The display shows one of the possible increments (1, 2, 5, 10, or 20 %), e.g. [10% Incr]. Select the desired value and press **CAL**.
4. Prepare a standard solution of the measured species. Its value should be 5 to 100 times higher than the expected sample concentration. The display shows one of the possible standards e.g. [Std 1.-5]. Select the desired value and press **CAL**.
5. After rinsing the electrodes with distilled water, immerse them in e.g. 100 ml sample solution. The display will show the electrode potential e.g. [E=178.2]. When readings are stable press **CAL**.
6. Add a small increment corresponding to the chosen mode (e.g. 5 ml). The display will show the difference in electrode potential e.g. [dE= 34.4]. When readings are stable press **CAL**. The instrument calculates the result and shows the concentration of the sample in the same unit as the standard used for the addition.
7. Calculate the real concentration by multiplying the result with the stoichiometric ratio (n). For a next measurement press **CAL** and the same procedure starts again from point 5). Press **MODE** to abort.

**Sample addition:**

*The analate addition method is used when the samples vary widely in temperature or have a 5 to 100 times higher concentration than the available standard. The most suitable increment can be found from: increment percentage < 100 x standard value / expected sample value.*

1. Press **MODE** until the display shows [Dir.CONC1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Anl.Add1]. Press **MODE** to proceed.
2. The display shows one of the possible increments (1, 2, 5, 10, or 20 %), e.g. [10% Incr]. Select the desired value and press **CAL**.
3. Prepare a standard solution of the measured species. Its value should be 5 to 100 times lower than the expected sample concentration. The display shows one of the possible standards e.g. [Std 3.-6]. Select the desired value and press **CAL**.
4. After rinsing the electrodes with distilled water, immerse them in e.g. 100 ml standard solution. The display will show the electrode potential e.g. [E= 281.5]. When readings are stable press **CAL**.
5. Add a small increment corresponding to the chosen mode (e.g. 20 ml). The display will show the difference in electrode potential e.g. [dE=23.9]. When readings are stable press **CAL**. The instrument calculates the result and shows the concentration of the sample in the same unit as the standard used for the addition.
6. For a next measurement press **CAL** and the same procedure starts again from point 4). Press **MODE** to abort.

**Sample subtraction:**

*The analate subtraction method is used to measure species for which stable standards or ion selective electrodes do not exist. The procedure is similar to the analate addition method. The most suitable increment can be found from: increment percentage < 100 x standard value / expected sample value.*

1. Calculate the stoichiometric ratio (n) of the reaction:  $xS + yR \rightarrow S_xR_y$   $n=x/y$  (S = sample, R = reagent).
2. Press **MODE** until the display shows [Dir.CONC1], [Std.Add1], [Std.Sub1], [Anl.Add1] or [Anl.Sub1] and choose [Anl.Sub1]. Press **MODE** to proceed.
3. The display shows one of the possible increments (1, 2, 5, 10, or 20 %), e.g. [10% Incr]. Select the desired value and press **CAL**.
4. Prepare a standard solution of the measured species. Its value should be 5 to 100 times lower than the expected sample concentration. The display shows one of the possible standards e.g. [Std 1.-6]. Select the desired value and press **CAL**.
5. After rinsing the electrodes with distilled water, immerse them in e.g. 100 ml standard solution. The display will show the electrode potential e.g. [E= 129.3]. When readings are stable press **CAL**.
6. Add a small increment corresponding to the chosen mode (e.g. 1 ml). The display will show the difference in electrode potential e.g. [dE= 41.7]. When readings are stable press **CAL**. The instrument calculates the result and shows the concentration of the sample in the same unit as the standard used for the addition.
7. Calculate the real concentration by multiplying the result with the stoichiometric ratio (n). For a next measurement press **CAL** and the same procedure starts again from point 5). Press **MODE** to abort.

**Maintenance:**

After prolonged use of the electrode the internal electrodes may become deactivated by undesired deposits. An indication is that the electrode no longer responds correctly to calibration. In this case:

1. Carefully screw off the measuring head.
  2. Cautiously remove the precipitated deposits with the aid of filter paper. **Do not use grinding paper or a glass fibre!**
  3. Rinse several times with distilled water and shake carefully off the water drops.
  4. Take the measuring head, fill it with new electrolyte solution and very slowly screw it onto the electrode while holding it vertically. **Take care that no air bubbles are present in the measuring head!**
  5. **Let the electrode rest for a few hours!** The electrode is now ready for use again.
- *Replace the membrane only when damaged, not when calibration is no longer possible!*

**Measurement:**

1. Select the mg/l (mg/l) or % range. The display will immediately show the measured oxygen according to the previous calibration. Should you want to recalibrate, press **CAL**.
  2. The display shows [SAL= MAn] (manual salinity correction) or [SAL= Aut] (automatic salinity correction). Choose the desired mode and press **CAL**. If you choose [SAL= MAn] the display will show, e.g. [SAL=21]. Leave this value to zero unless you are going to measure in heavily salted solutions such as e.g. sea-water (35). Pre-set the proper value and press **CAL**.
  3. The display shows the pressure correction, e.g. [P=1023hP]. Pre-set the actual air pressure (hP) and press **CAL**.
  4. The electrode exposed to the air reaches an equilibrium corresponding to the partial pressure of oxygen and thus to saturation in water at the given temperature. The instrument shows the measured saturation, e.g. [SA<sub>t</sub>=8.3] and will calibrate automatically when readings are stable ([=] stops blinking).
  5. After rinsing the electrodes with distilled water, immerse them in the samples and read the display. The advection rate must be at least 10 cm/s (**stir!**).
  6. Rinse the electrode always after use and store it in distilled water.
- *Press RES to change the resolution from 0.1 to 0.01 mg/l or from 1 to 0.1 %.*

**Relative O<sub>2</sub> measurement:**

1. While in the O<sub>2</sub> range, press **▲**. At that moment the instrument stores the actual value and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
2. To cancel relative readings, press **MODE** and you will be in the normal measuring mode again.

1. Press **SET** until the display shows [diAGn].
2. Sequentially press **CAL** to show the calibration details of all connected electrodes, e.g:  
[94.6%] (slope)  
[6.89 pHo] (ISO-potential)  
[cc=1.034] (cell constant)  
[Cur=67.9] (saturation-current (µA))  
[Sat=8.73] (saturation-point (mg/l))
3. Press **SET** to return or press **PRINT** each time to print a complete calibration report of the selected channel, e.g.:

## Multi-channel Analyser C835

```
-----  
Channel           : pH1  
Date              : 21/09/95  
Time              : 10:06:01
```

## SETTINGS

```
-----  
Identification No : 000  
Software Version  : 1.0  
User Code         : off
```

```
ATC                (Pt1000) : on  
Calibr. Temp.      (C)       : 23.4  
Ref. Temp.         (C)       : 25.0
```

## CALIBRATION

```
-----  
Date               : 21/09/95  
Time               : 10:05:26
```

```
Buffer 1.68 / 4.01  
Slope              (%)       : 98.1  
Zero Point         (pH)      : 6.85
```

```
Buffer 4.01 / 6.87  
Slope              (%)       : 99.0  
Zero Point         (pH)      : 6.89
```

```
Buffer 6.87 / 9.18  
Slope              (%)       : 97.3  
Zero Point         (pH)      : 6.87
```

```
Buffer 9.18 / 12.45  
Slope              (%)       : 96.2  
Zero Point         (pH)      : 6.81
```

```
Average Slope     (%)       : 97.7  
Average pHo       (pH)      : 6.86  
Average Resp.     (s)       : 14  
Average Temp.     (C)       : 21.9
```

## STATISTICS

```
-----  
Average Slope     (%)       : +1.8  
Average pHo       (pH)      : +0.31  
Average Resp.     (s)       : -1
```

**Pre-set and start the data-logging:**

1. Press **SET** until the display shows [SEt LoG] and press **CAL**.
2. Select [Log dAtA] and press **CAL**.
3. Choose between [LOG PH], [LOG MV], [LOG MS],... etc. Select the desired mode and press **CAL**.
4. The display shows e.g. [dt=100]. Select the desired time interval (0...9999 s) between the data-logging and press **CAL**.
5. This step is **skipped** unless manual data-logging is chosen. Pre-set the time-interval (dt) to zero and press **CAL** (point 6 is skipped).
6. The display shows e.g. [n=40]. Select the desired number of values (1...500) to be data-logged and press **CAL**.
7. The display shows [StArt ?]. Press **CAL** to start the data-logging according the previous settings (a blinking [L] appears) or press **MODE** to return. After the last value has been data-logged the display will show [End LOG]. When manual data-logging has been selected, press **PRINT** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [LOG 7]. Press **MODE** to return.

**Continue with a next series of measurements, without erasing the previous data-logging:**

1. Press **SET** until the display shows [SEt LoG] and press **CAL**.
2. Select [Cnt LoG] provided the memory has enough capacity left! Press **CAL** to continue with a next series of measurements, without erasing the previous data-logging.
3. Press **SET** to return.

**View the stored values on the display:**

1. Press **SET** until the display shows [SEt LoG] and press **CAL**.
2. Select [LSt dAtA] and press **CAL**.
3. The display shows the first data-logged value in memory (a blinking [n] appears). Press  $\blacktriangle$  or  $\blacktriangledown$  to view all other values.
4. While pressing **HOLD** the corresponding date is shown.
5. While pressing **RES** the corresponding time shown.
6. Press **PRINT** to send the recalled data to a printer.
7. Press **SET** to return.

**Send the stored values to a printer or computer:**

1. Press **SET** until the display shows [SEt LoG] and press **CAL**.
2. Select [Prt dAtA] and press **CAL** to print the data.
3. The display shows [PrintinG] while all data-logged values, including corresponding dates and times, are sent to the RS232 output in the following format, compatible with EXCEL™, LOTUS 123™,....The columns are separated by a TAB character:

Identification No: 037

#	VALUE	UNIT	C	CH	H	D
0001	7.34	pH	23.7	2	13:19:57	29/01/95
0002	7.21	pH	23.7	2	13:20:57	29/01/95
0003	7.08	pH	23.8	2	13:21:57	29/01/95
0004	6.71	pH	23.8	2	13:22:57	29/01/95

**Interfacing with a computer:**

Send a single character to the instrument and it will execute the corresponding command as follows:

- 1 = press **MODE**.
- 2 = press **CAL**.
- 3 = press **▲**.
- 4 = press **▼**.
- 5 = press **SET**.
- 6 = press **HOLD**.
- 7 = press **RES**.
- 8 = press **PRINT**.
- ? = send display to computer.
- + = keyboard on (manual operation remains possible).
- = keyboard off (manual operation disabled).
- V = start the input of a value (only valid in certain routines).

Follow exactly the same measuring or calibration procedures, as described in this manual, to include any desired command in your computer data-acquisition program. This example continuously prints the measured values on the computer screen:

```

10  ON ERROR GOTO 70           'makes sure no error occurs
20  OPEN "COM1:2400,N,8,2" AS #1 'prepares computer input
30  PRINT #1,"8"              'PRINT button pressed
40  INPUT #1,V$               'measured values transmitted
50  PRINT V$                  'received values printed
60  GOTO 30                   'repeat procedure
70  RUN

```

**Input of a value:**

The direct input of values ("V" command) is only possible when the instrument is running a routine where also a manual input of these values could be made. If not, it will ignore any "V" command from the computer!

**Examples of allowed routines are:**

- temperature input (e.g. manual temperature compensation).
- control parameter input such as Low, High, ...
- time settings.
- input of manual buffer values.

**Examples of NOT allowed routines are:**

- during measurements.
- when values have to be selected rather than changed (e.g. buffer values in memory).
- while an electrode is being calibrated.

A 16-bit value (2's complement) should be transmitted in the following sequence:

```

1st character = "V" (start a value input)
2nd character = highest byte of the value in ASCII
3rd character = lowest byte of the value in ASCII
4th character = checksum of 2nd and 3rd character in ASCII
5th character = LF, linefeed (ASCII-10)

```

When a correct checksum has been received, the instrument will send a confirmation character "!" to the computer. If not, a question mark "?" is sent. Both are eventually preceded by the identification number. This example makes the computer to enter the value "1000" on instrument #7, in the sophisticated mode:

```

1000 = (H-byte 3)(x 256) + (L-byte 232)
10  OPEN "COM1:2400,N,8,2" AS #1   'prepares communication port
20  PRINT #1,"V";                  'start a value input
30  PRINT #1,CHR$(3);              'highest byte of value
40  PRINT #1,CHR$(232);            'lowest byte of value
50  PRINT #1,CHR$(235);            'checksum of both bytes
60  PRINT #1,CHR$(10);             'LF, closes command

```

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If you need assistance with a technical problem, make sure to fill the **complete** checklist (all 23 points!) and fax a copy to Consort: +32/14/429179.

1. purchase date:

2. purchased from:

2. indicate model:

C831 - C832 - C833 - C835 - C838

4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code:

10. city:

11. telephone:

12. telefax:

13. e-mail:

14. type of connected electrode:

**Perform the following steps:**

15. store all pH electrodes in a 3...4 KCl solution overnight.

yes?

16. reset the meter (switch on while holding MODE pressed).

yes?

17. short-circuit the pH input, select the mV mode and read display.

mV:

18. select the °C mode, measure temperature of solutions and read the display.

°C:

19. select the mV mode (do not calibrate!).

yes?

20. connect the pH electrode, rinse, dip in fresh buffer 7 (stir!) and read display.

mV:

21. rinse the pH electrode, dip in fresh buffer 4 (stir!) and read display.

mV:

22. connect a new pH electrode, rinse, dip in fresh buffer 7 (stir!) and read display.

mV:

23. rinse this pH electrode, dip in fresh buffer 4 (stir!) and read display.

mV:

If you need assistance with a technical problem, make sure to fill the complete checklist (all 23 points!) and fax a copy to Consort: +32/14/429179.

1. purchase date:  2. purchased from:

2. indicate model:  4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code:  10. city:

11. telephone:  12. telefax:

13. e-mail:

14. type of connected electrode:

Perform the following steps:

- |  |      |                          |
|--|------|--------------------------|
| 15. store all conductivity cells in distilled water with some detergent overnight.       | yes? | <input type="checkbox"/> |
| 16. reset the meter (switch on while holding MODE pressed).                              | yes? | <input type="checkbox"/> |
| 17. leave the EC input open, select the conductivity mode and read display.              | µS:  | <input type="checkbox"/> |
| 18. connect a 1000 ohm resistance to the conductivity input and read display.            | µS:  | <input type="checkbox"/> |
| 19. select the °C mode, measure temperature of solutions and read the display.           | °C:  | <input type="checkbox"/> |
| 20. rinse the conductivity cell with a 0.01 M KCl calibration solution (1413 µS).        | yes? | <input type="checkbox"/> |
| 21. connect the conductivity cell, dip in fresh cal. solution (stir!) and read display.  | µS:  | <input type="checkbox"/> |
| 22. rinse a new conductivity cell with a 0.01 M KCl calibration solution (1413 µS).      | yes? | <input type="checkbox"/> |
| 23. connect this conductivity cell, dip in fresh cal. solution (stir!) and read display. | µS:  | <input type="checkbox"/> |

If you need assistance with a technical problem, make sure to fill the **complete** checklist (all 23 points!) and fax a copy to Consort: +32/14/429179.

1. purchase date:  2. purchased from:

2. indicate model:  4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code:  10. city:

11. telephone:  12. telefax:

13. e-mail:

14. type of connected electrode:

**Perform the following steps:**

- |  |      |                          |
|--|------|--------------------------|
| 15. prepare the ion electrode (ISE) and solutions according to the manual.               | yes? | <input type="checkbox"/> |
| 16. reset the meter (switch on while holding MODE pressed).                              | yes? | <input type="checkbox"/> |
| 17. short-circuit the pH input, select the mV mode and read display.                     | mV:  | <input type="text"/>     |
| 18. select the °C mode, measure temperature of solutions and read the display.           | °C:  | <input type="text"/>     |
| 19. connect the ISE, select the mV mode (do <b>not</b> calibrate!).                      | yes? | <input type="checkbox"/> |
| 20. rinse the ISE, dip in a fresh 1 ppm cal. solution + ISA (stir!) and read display.    | mV:  | <input type="text"/>     |
| 21. rinse the ISE, dip in a fresh 10 ppm cal. solution + ISA (stir!) and read display.   | mV:  | <input type="text"/>     |
| 22. rinse the ISE, dip in a fresh 100 ppm cal. solution + ISA (stir!) and read display.  | mV:  | <input type="text"/>     |
| 23. rinse the ISE, dip in a fresh 1000 ppm cal. solution + ISA (stir!) and read display. | mV:  | <input type="text"/>     |

If you need assistance with a technical problem, make sure to fill the complete checklist (all 20 points!) and fax a copy to Consort: +32/14/429179.

1. purchase date:

2. purchased from:

2. indicate model:

C831 - C832 - C833 - C835 - C838

4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code:

10. city:

11. telephone:

12. telefax:

13. e-mail:

14. type of connected electrode:

Perform the following steps:

15. Carefully screw off the measuring head of the oxygen electrode. Cautiously remove the precipitated deposits with the aid of filter paper. Do not use grinding paper or a glass fibre! Rinse several times with distilled water and shake carefully off the water drops. Take the measuring head, fill it with new electrolyte solution and very slowly screw it onto the electrode while holding it vertically. Let the electrode rest for a few hours!      yes?

16. reset the meter (switch on while holding MODE pressed).      yes?

17. short-circuit the O<sub>2</sub> input, select the O<sub>2</sub> mode and read display.      mg/l:

18. select the °C mode, measure air temperature and read the display.      °C:

19. select the O<sub>2</sub> mode (do not calibrate!).      yes?

20. rinse the O<sub>2</sub> electrode, hold in the air and read display.      mg/l:

We would like to know your opinion, provide us with your remarks and fax a copy of this page to Consort: +32/14/429179.

1. purchase date:

2. purchased from:

2. indicate model:

4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code:

10. city:

11. telephone:

12. telefax:

13. e-mail:

- Are you generally satisfied with the performance of the instrument ?
- Are you satisfied with the service of the company you purchased from ?
- What is the degree of satisfaction with:
  - the design of the instrument ?
  - the operating instructions ?
  - the supplied accessories ?
  - the delivery ?
  - the price/quality ratio ?
- What are your suggestions:

		
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....

.....