CONSORT

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

E122

E132

E143

QUALITYMASTERS

CERTIFICATE OF REGISTRATION

This is to certify that:

Consort NV

Parklaan 36
B-2300 Turnhout

Has been assessed by QualityMasters in respect of its Quality Management System and found to comply with:

NEN-EN-ISO 9001 : 2000

Approval is hereby granted for registration providing the rules and conditions relating to certification are observed at all times.

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Lack of fulfillment of conditions as set forth in the certification regulations may render this certificate invalid. The use of the accreditation mark indicates accreditation in respect to the activities covered by the accreditation number EU9706010F.

Q

M

REGISTRATION

CERTIFICATE OF

Authorized signatory

A. LEUWATY
QualityMasters B.V.

This certificate remains the property of QualityMasters B.V.

WARRANTY CERTIFICATE

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory.

CONSORT will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse.

A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

DECLARATION OF CONFORMITY

We declare under our sole responsibility that the product

Electrophoresis Power Supply

content of the type numbers

E122, E132, E143

to which this declaration relates is in conformity

with the following standards

EN61010

LOW VOLTAGE DIRECTIVE 73/23/EEG

EN50081-1

EN50082-1

EN60555-2

EMC DIRECTIVE 89/336/EEG

Turnhout, September 8, 1998

A handwritten signature in black ink, appearing to be 'J. G. G.', is written over a horizontal line.

on behalf of **CONSORT nv**

This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 4° and 40° C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

Connect the instrument only to an earthed power line. The required power source is indicated on the label at the back of the instrument. Do not cut and splice the power cord. When removing the power cord from the wall outlet, be sure to unplug by holding the plug attachment and not by pulling the cord. Do not hold the plug by wet hand.

Manufacturer

CONSORT nv

Parklaan 36
B2300 Turnhout
Belgium

Tel (+32)(14)41 12 79
Fax (+32)(14)42 91 79
E-mail: info@consort.be

Warranty

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

CONSORT is not liable for consequential damages arising out of the use or handling of its products.

Servicing

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply lead and **NOT** to send the following items unless they are suspect:

Manual

Cables

Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

Ranges	E122: 0...150 V, max. 150 mA E132: 0...300 V, max. 150 mA E143: 0...400 V, max. 250 mA
Outputs	3 pair of 4 mm sockets
Ambient temp.	4...40°C
Relative humidity	0...90 % (non-condensing!)
Power supply	210-250 V~ (105-125 V~), 50/60 Hz
Cabinet	rugged plastic cabinet
Dimensions	15 x 20 x 12 cm
Weight	2 kg

Very important

This electrophoresis power supply is a high technology instrument available in several versions. As it is capable of giving dangerous voltage levels by which high power is involved, we suggest that you take a few moments to read this manual thoroughly. Although this instrument is equipped with all necessary safety features against abuse and other accidental failures, caution should be exercised when working with high voltage equipment. Therefore, **avoid to touch the outlets with any conducting object** and make sure there is a second person present for your safety in case of any severe electric shock. Never touch any part of the assembly (power supply, leads or tank) before having switched off. **Never manipulate with wet hands. Do not ground any of the outputs or the buffer in the tank.** Connect the outlets only to an insulated electrophoresis tank with safety cover. Never make any other connections, such as e.g. putting several power supplies in series or in parallel. In order to prevent electric shock, never open the back plate nor remove the cover. Do not expose the unit to rain or any other liquid. **Do not spill liquid or insert metal objects inside the unit.** Take care so that the power supply is not dropped to avoid damaging the cabinet which defeats safeguards or injuring yourself. If the unit has been dropped or the cabinet has been damaged, unplug it and have it checked by an authorised service technician to restore the safeguards. The fact that the unit operates satisfactorily does not imply that the unit is properly earthed or that it is completely safe. If in any doubt about the effective earthing of the unit, contact a qualified electrician. **Never block the ventilation holes** or place the unit in any enclosure unless proper ventilation is provided. Never place the unit near or over a radiator, heat register or stove. Avoid locations where the instrument is exposed directly to the sun light.

Operation

1. Verify if the instrument is switched off.
2. Connect the electrophoresis tank to the instrument.
3. Turn the voltage adjuster to zero.
4. Switch the instrument on.
5. Turn the voltage adjuster slowly to the desired voltage.

When the alarm indicator illuminates, the maximum current is exceeded and the instrument will automatically limit the voltage to a safe value.

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If you need assistance with a technical problem, make sure to fill the **complete** checklist (26 points!) and fax a copy to Consort: +32/14/429179.

1. purchase date: 2. purchased from:

2. indicate model: 4. serial no:

5. organisation:

6. name:

7. division:

8. address:

9. postal code: 10. city:

11. telephone: 12. telefax:

13. e-mail:

14. type of connected electrophoresis unit:

Perform the following steps:

15. connect the electrophoresis unit (or a dummy resistance)	yes?	<input type="checkbox"/>
16. turn the voltage button to the desired value.	V:	<input type="text"/>
17. read the output current.	mA:	<input type="text"/>
18. does the alarm sign illuminate.	yes/no?	<input type="checkbox"/>
19. leave ALL outputs open (no load connected).	yes?	<input type="checkbox"/>
20. turn the voltage button to the desired value.	V:	<input type="text"/>
21. connect a voltmeter between the red & black outlets and measure the voltage.	V:	<input type="text"/>