# CONSORT

---

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

**K611**

---

January 1997

## **Table of contents**

Introduction .....	1
Manufacturer .....	1
Warranty .....	1
Servicing .....	1
Specifications .....	2
Keyboard .....	2
Error Codes .....	3
AC Adapter .....	3
<b>SET-UP .....</b>	<b>3</b>
Alarm .....	3
Beeper .....	4
Printer .....	4
Identification .....	4
Code .....	4
Temperature compensation .....	4
<b>DATA ACQUISITION .....</b>	<b>5</b>
Logging data .....	5
Adding data .....	5
Recalling data .....	6
Printing data .....	6
Inputs .....	6
<b>CONDUCTIVITY MEASUREMENT .....</b>	<b>6</b>
Maintenance .....	6
Calibration .....	6
Conductivity measurement .....	7
<b>TEMPERATURE MEASUREMENT .....</b>	<b>7</b>
Calibration .....	7
Measurement .....	7
<b>DATA COMMUNICATION .....</b>	<b>8</b>
Recorder output .....	8
GLP-protocol .....	8
Digital output .....	9
Printer connection .....	9
Computer connection .....	9
<b>TROUBLE SHOOTING .....</b>	<b>10</b>
Simple test .....	10
Calibrate .....	10
Reset .....	10

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This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 4° and 40°C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

If it is equipped with rechargeable batteries, the instrument should be connected permanently to the mains in order to keep the batteries in good condition. Do not store the instrument for longer than 1 month without recharging the batteries. The batteries can be replaced by opening the bottom compartment of the cabinet.

**CONSORT nv**  
Parklaan 36  
B2300 Turnhout  
Belgium

Tel (++32)(14)41 12 79  
Fax (++32)(14)42 91 79

## ***Introduction***

## ***Manufacturer***

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

**CONSORT** is not liable for consequential damages arising out of the use or handling of its products.

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply and **NOT** to send the following items unless they are suspect:

*Manual*  
*Cables*  
*Accessories*

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

## ***Warranty***

## ***Servicing***

## Specifications

<b>Conductivity</b>	0...200 mS/cm
<b>Resistivity</b>	0...20 MΩ.cm
<b>Temperature</b>	0...100°C
<b>Resolution</b>	0.1°C, 0.01 μS/cm, 10 Ω.cm
<b>Inputs</b>	BNC input for conductivity cell BANANA input for a Pt1000 temperature probe
<b>Temp. Comp.</b>	automatic with Pt1000 or manual
<b>Digital output</b>	programmable RS232, 2400 b/s, for bi-directional communication with a computer or printer
<b>Analogue output</b>	0...2 V, ca 8 kΩ, 8 bit D/A
<b>Display</b>	8 digit/character, 12 mm LCD display, shows readings, °C & messages
<b>Keys</b>	10 tactile membrane keys
<b>Ambient temp.</b>	4...40°C
<b>Rel. Humidity</b>	0...90 % (non-condensing!)
<b>Power supply</b>	210-250 V~ (* 105-125 V~), 50/60 Hz, max. 2 VA, & 4 NiCd batteries, size AA
<b>Cabinet</b>	IP65 cabinet
<b>Dimensions</b>	252 x 121 x 50 mm
<b>Weight</b>	600 g

Each time the instrument is switched on, it will check briefly its internal calibrations while showing all functions of the display and the software-version. An error message appears when something is wrong.

<b>MODE</b>	= Selects all modes or escapes from error traps, calibration procedures, etc..by returning to the original mode.
<b>CAL</b>	= Starts or proceeds a calibration or a function.
<b>▲/▼</b>	= Button for entering a value or for selecting a function.
<b>°C</b>	= Reads temperature, when a Pt1000 is present, or manual temperature compensation.
<b>HOLD</b>	= Holds display when measuring.
<b>RES</b>	= Changes resolution/units of the display.
<b>PRINT</b>	= Displayed value is printed through the RS232 output.
<b>ON/OFF</b>	= Switches the instrument on or off.

## Keyboard

## **Error codes**

- [Or. EC] = Overrange Conductivity (e.g. short circuited electrode cable).
- [Or. OhM] = Overrange Resistivity (e.g. interrupted electrode cable).
- [Or. °C] = Overrange °C (measured degrees out of range).
- [no dAtA] = Data-logging memory empty.
- [not.CAL] = Calibration procedure interrupted (the previous calibration remains valid!).
- [Err.cc] = Cell constant not within 70...130% of the nominal value.
- [Err.CAL] = Calibration error (check electrodes and standard solutions).
- [Err.Stb] = Electrode not stable.
- [Err.Prt] = Communication error with the attached printer.
- [Err.bAt] = Battery charger defective (disconnect the AC-adapter immediately from the mains, remove the batteries and have your instrument serviced!).
- [Err.COM] = General communication error.
- [Err.MEM] = Fatal memory error (all factory settings are cleared. Have your instrument serviced!).
- [ILLEGAL] = An unauthorised person has failed to repair the unit.  
**Warranty is elapsed from now on!**

Remove the protective silicone stopper from the REF socket and connect the jack of a suitable AC adapter (model **A4070** for 230 V~ or model **A4071** for 115 V~) to the socket for recharging the NiCd batteries. **Avoid to use an AC adapter without NiCd batteries being inserted in the instrument!** For field work, remove the adapter and reinsert the protective silicone stopper into the socket.

## **AC Adapter**

# SET-UP

*An audible alert can be given when the displayed values stray outside a programmable high- and low-level.*

## **Alarm**

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [ALrM OFF], [ALrM EC], or [ALrM tdS]. Press **CAL**. Select [ALrM OFF] if no alarm is required.
3. The display shows e.g. [L=23.0MS] while the units are blinking. Select the desired range with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**.
4. The display shows e.g. [L=23.0MS] while [=] is blinking. Select the desired minimum level with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Each time the measurements are below this level, the beeper will sound an alarm.
5. The display shows e.g. [H=57MS] while the units are blinking. Select the desired range with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**.
6. The display shows e.g. [H=57MS] while [=] is blinking. Select the desired maximum level with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Each time the measurements are above this level, the beeper will sound an alarm.
7. Press **CAL** to continue or press **MODE** to cancel.

## **Beeper**

*The built-in beeper can be programmed to sound each time a key-path is pressed.*

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [bEEP on] or [bEEP OFF]. Decide whether or not the beeper should sound each time the keyboard is touched.
3. Press **CAL** to continue or press **MODE** to cancel.

## **Printer**

*The measured values can be printed automatically at timed intervals or manually on operator command.*

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Pt=120]. Select the desired time interval (0...9999 s) between the transmitted data to a printer, with  $\blacktriangle$  or  $\blacktriangledown$ . Pre-set to zero if no automatic printing is required.
3. Press **CAL** to continue or press **MODE** to cancel.

## **Identification number**

*The instrument can be identified by e.g. a computer when a specific number is allocated to it.*

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Id.no=5]. Select the desired identification number (0...999) for your meter with  $\blacktriangle$  or  $\blacktriangledown$ . Pre-set to zero if this number should not be mentioned while printing.
3. Press **CAL** to continue or press **MODE** to cancel.

## **Secret code**

*A private code can be programmed to avoid undesired access to the instrument.*

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows [Code on], [Code OFF], or [Code CAL]. Decide, with  $\blacktriangle$  or  $\blacktriangledown$ , whether the access to all future calibrations should be by entering a code (on) or not (off). Press **CAL** to continue or press **MODE** to cancel. Select [Code CAL] to enter your personal code.
3. The display shows [E=-----] while [=] is blinking. Enter your secret sequence of 5 keys to proceed. Press **CAL** to continue or press **MODE** to cancel.

## **Temperature compensation**

*Permits to install the reference temperature.*

1. Select the set-up procedure by pressing **MODE**.
2. The display shows [rEF=25°C] or [rEF=20°C]. Select the temperature to which all future conductivity measurements will be referred with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL** to continue or **MODE** to return.

# DATA-ACQUISITION

*Up to 100 values can be stored in the non-volatile memory of the instrument.*

- [no LOG]** : No data-logging.
- [SEt LOG]** : Permits to pre-set the data-logging parameters.
- [LSt LOG]** : Permits to view the stored values on the display.
- [Prt LOG]** : Permits to send the stored values to a printer or computer.
- [Cnt LOG]** : Permits to continue with a next series of measurements, without erasing the previous data-logging.

## **Set log: Procedure to pre-set the data-logging parameters.**

## ***Logging data***

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [SEt LOG] with  $\blacktriangle$  or  $\blacktriangledown$ . Select [no LOG] if no data-logging is required and press **MODE** to return to the original mode.
3. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [LOG EC], [LOG OhM] or [LOG tdS]. Select the desired mode and press **CAL**.
4. The display shows e.g. [dt=100]. Select the desired time interval (0...9999 s) between the data-logging, with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Pre-set to zero for manual data-logging, each time **PRINT** is pressed (the next points 5 to 7 are skipped and press **CAL**).
5. The display shows e.g. [n=40]. Select the desired number of values (1...100) to be data-logged with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**.
6. The display shows e.g. [d=31.12.91] while the year is blinking. Program the desired starting date with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Program similarly the month and the day.
7. The display shows e.g. [t=23.59] while the minutes are blinking. Program the desired starting time with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Program similarly the hours.
8. The display shows [StArt ?]. Press **CAL** to start the data-logging according the previous settings (a blinking [L] appears) or press **MODE** to cancel. After the last value has been catalogued the display will show [End LOG]. When manual data-logging has been selected, press **PRINT** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [LOG 7]. Press **MODE** to cancel.

## **[Cnt LOG]: Permits to continue with a next series of measurements, without erasing the previous data-logging**

## ***Adding data***

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Cnt LOG] with  $\blacktriangle$  or  $\blacktriangledown$  to continue with a next series of measurements, without erasing the previous data-logging, provided the memory has enough capacity left and press **CAL**.

**[Lst LOG]: permits to view the stored values on the display.**

## **Recalling data**

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Lst LOG] with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**
3. The display shows the first data-logged value in memory (a blinking [n] appears). Press  $\blacktriangle$  or  $\blacktriangledown$  to view all other values.
4. While pressing **HOLD** the corresponding date is shown (not relevant when a manual or continued data-logging has been performed).
5. While pressing **RES** the corresponding time shown (or: series + log-number, e.g. [2-017], when a continued data-logging has been performed).
6. While pressing **°C** the corresponding temperature is shown.
7. Press **PRINT** to send the recalled data to a printer.
8. Press **MODE** to return to cancel.

**[Prt LOG]: Permits to send the stored values to a printer or computer.**

## **Printing data**

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [Prt LOG] with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**
3. The display shows [PrintinG] while all data-logged values, including corresponding dates and times, are sent to the RS232 output in the following format, compatible with EXCEL™, LOTUS 123™,... The columns are separated by a TAB character.

Log#	VALUE	UNIT	C	H	D
001	16.49E-03	S/cm	23.7	13:19:57	29/01/95
002	15.05E-03	S/cm	23.7	13:20:57	29/01/95
003	13.39E-03	S/cm	23.8	13:21:57	29/01/95

The measuring cell should be connected to the coaxial INPUT connector. Automatic temperature compensation and temperature measurements are possible by plugging a Pt1000 thermocompensator into the °C terminals. You can also use a combination conductivity electrode with built in Pt1000. Its banana plug should be inserted in the upper °C terminal. Without compensator, the manual temperature compensation is automatically switched on.

## **Inputs**

# **CONDUCTIVITY MEASUREMENT**

**A conductivity cell is active and stable only after wetting!** For this purpose it must be immersed for **at least one hour** in distilled water. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

## **Maintenance**

A polluted cell may be cleaned with a soft detergent or diluted nitric acid. Greasy substances may be removed with acetone or alcohol (never do this with plastic electrodes!).

*The cell constant depends on the physical dimensions of the platinum plates. They are coated with a spongy black platinum deposit. This increases greatly the effective surface and reduces polarising effects. Since the coating of platinum black is quite resistant to chemical action, the cell constant remains fairly constant during the electrode life time.*

## **Calibration**

1. Select the conductivity range by pressing **MODE** (the display shows [ $\mu$ S] or [MS]) and press **CAL**.
2. The display will show the cell constant (e.g. [cc=1.073]). Press **CAL** to start the automatic determination procedure.
3. The display shows one of the KCl standard solutions in memory [St=0.01M] (1413  $\mu$ S/cm), [St=0.1M] (12.88 mS/cm) or [St=1M] (111.8 mS/cm). Select the proper standard with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**.
4. After rinsing the cell several times with the standard solution, immerse it the standard solution. The solution temperature is not so critical but it should lie between 0°C and 30°C. If the standard has a different temperature press  $^{\circ}$ **C** here (not necessary when using a Pt1000). Press **CAL** to continue.
5. The display will show the computed cell constant e.g. [cc1=1.045] and will calibrate automatically when readings are stable ([=] stops blinking).

1. Select the conductivity range by pressing **MODE** (the display shows [ $\mu$ S] or [MS]).
2. Rinse the electrodes with the sample and immerse them in it.
3. Read the display (remember: compensated to the reference temperature!). The autoranging feature gives you the largest possible number of digits + their unit.
4. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

## ***Conductivity measurement***

- \* Disconnect the AC-adapter to increase accuracy in the lower conductivity range!
- \* See page 41 for the standards in memory.
- \* A blinking decimal point warns you for unstable measurements. Wait to read the display!
- \* Press **RES** to switch between ppm (salinity), S/cm (conductivity) or  $\Omega$ .cm (resistivity).
- \* When starting a titration, press  $\blacktriangledown$  to lock the actual range and avoid cross-over errors due to differences in the measuring frequency. Press **MODE** to return to the normal measuring mode.
- \* Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.

# **TEMPERATURE MEASUREMENT**

1. Press **CAL** to calibrate the  $^{\circ}$ C scale with any reference. Immerse the Pt1000 in a solution of known temperature.
2. The display shows e.g. [ $^{\circ}$ C=22.3]. Calibrate to the proper value with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL** again or press **RES** to reset the absolute readings.

## ***Calibration***

1. Read temperature by pressing  $^{\circ}$ **C**.
2. Without Pt1000, adjust the manual temperature compensation with  $\blacktriangle$  or  $\blacktriangledown$  and proceed by pressing **MODE**.

## ***Measurement***

# DATA COMMUNICATION

A recorder can be connected to the red (+) and the black (-) terminals. The 0 to 2 V output range corresponds with 0 to 2000 digits. Use only laboratory recorders with a high input impedance!

## **Recorder output**

1. Select the conductivity range by pressing **MODE**.
2. Press **CAL** to display the cell constant, e.g. [cc=1.073].
3. Press **PRINT** to print a complete calibration report, e.g.:
4. Press **MODE** to return to the normal measuring mode.

## **GLP-protocol**

```
CONSORT Conductometer K611
-----

SETTINGS
-----
Identification No      : 004
Software Version      : 3.0
User Code              : on
Titration Lock        : off

ATC      (Pt1000)      : yes
ATC Cal. Temp. (C)    : 30.0

Alarm EC              : on
Alarm High (S/cm)     : 85.7E-03
Alarm Low  (S/cm)     : 30.5E-03

Ref. Temp.            : 20

Batt.Capacity (%)     : 87

CALIBRATION
-----
Ref. Temp. (C)        : 25
Standard (S/cm)       : 1413E-06
Cell Const (cm-1)     : 1.073
Response Time (s)     : 13

STATISTICS
-----
Cell Const (cm-1)     : -0.026
Average Resp. (s)     : +1
```

## Digital output

A standard RS232 output terminal (DP9) is provided for interfacing the instrument with a printer or computer. Data is sent in the ASCII code at a BAUD rate of 2400 b/s (8 bit, no parity, 1 start and 2 stopbits, automatic line feed ON, data transfer protocol Xon/Xoff).

### Serial port pinout specifications:

pin 1 : RLSD, received line signal detector  
pin 2 : TxD, transmit data  
pin 3 : RxD, receive data  
pin 4 : not connected  
pin 5 : Gnd, signal ground  
pin 6 : DSR, data set ready  
pin 7 : not connected  
pin 8 : CTS, clear to send  
pin 9 : not connected

### Interfacing with a printer:

Press **PRINT** each time you want to print the measured values. The columns are separated by a TAB character. Run the set-up procedure (page for automatic printing at timed intervals, 'Printer' page 4, while you are measuring. The **PRINT** keypath does not function in this case. It is also disabled during an automatic data-logging., e.g.:

Log#	VALUE	UNIT	C	H	D
001	16.49E-03	S/cm	23.7	13:19:57	29/01/95
002	15.05E-03	S/cm	23.7	13:20:57	29/01/95
003	13.39E-03	S/cm	23.8	13:21:57	29/01/95

## Printer connection

### Using WINDOWS™ to collect data:

1. Connect the computer with the RS232 terminal, using the optional AK2310 cable (eventually with adapter AB3190 for connection with a 25 pole RS232 terminal).
2. Start **WINDOWS™** and select *Terminal* in the *Accessories Group*.
3. Choose *Receive Text File* in the *Transfer Menu* and create a file name (TXT format) for the data to receive.
4. Select the *Settings Menu* and make sure the following parameters are programmed:

Baud Rate : 2400  
Data bits : 8  
Stop bits : 2  
Parity : None  
Flow control : Xon/Xoff

5. Start the conductometer in the desired measuring mode for manual or automatic data collection (see 'Printer' page 4 for more details).
6. When finished, open the created file with a standard spreadsheet (e.g. **EXCEL™**) or wordprocessor (e.g. **WORD™**) for further processing.

## Computer connection

## Interfacing with a computer:

Send a single character to the instrument and it will execute the corresponding command as follows:

- 1 = press **MODE**.
- 2 = press **CAL**.
- 3 = press **▲**.
- 4 = press **▼**.
- 5 = press **SET**.
- 6 = press **HOLD**.
- 7 = press **RES**.
- 8 = press **PRINT**.
- ? = send display to computer.
- + = keyboard on (manual operation remains possible).
- = keyboard off (manual operation disabled).

Follow exactly the same measuring or calibration procedures, as described in this manual, to include any desired command in your computer data-acquisition program.

This example continuously prints the measured values (pH and temperature) on the computer screen:

10	ON ERROR GOTO 70	'makes sure no error occurs
20	OPEN "COM1:2400,N,8,2" AS #1	'prepares computer input
30	PRINT #1,"8"	' <b>PRINT</b> button pressed
40	INPUT #1,V\$	'measured values transmitted
50	PRINT V\$	'received values printed
60	GOTO 30	'repeat procedure
70	RUN	

# TROUBLE SHOOTING

*Try this procedure if you suspect the electronic performance of the instrument.*

1. Leave the conductivity input open.
2. Switch the instrument ON.
3. Select the conductivity range by pressing **MODE**. The display should show  $[0.00\mu\text{S}] \pm 0.01 \mu\text{S/cm}$ . Run the **Calibrate** procedure if erroneous values are read.

*This procedure permits to calibrate automatically all amplifiers of the instrument.*

1. Leave the conductivity input open.
2. Switch the instrument ON while holding **CAL** pressed.
3. The display shows [SELF CAL] for a while and returns to the measurements.

*This procedure resets the instrument to the original factory settings. All personal memorised data (including access code) will be erased!*

1. Switch the instrument ON while holding **MODE** pressed.
2. The display shows [rESET] for a while and returns to the measurements.

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