CONSORT

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

K612

January 1997

Table of contents

Introduction	1
Manufacturer	1
Warranty	1
Servicing	1
Specifications	2
Keyboard	3
Error Codes	3
AC Adapter	3
SET-UP	4
Beeper	4
Alarm	4
Printer	4
Identification	4
Code	5
Temperature compensation	5
DATA ACQUISITION	5
Logging data	5
Adding data	6
Recalling data	6
Printing data	6
Inputs	7
CONDUCTIVITY MEASUREMENT	7
Maintenance	7
Calibration	7
Temperature compensation	8
Capacitance compensation	8
Conductivity measurement	9
Salinity measurement	9
Relative measurement	9
TEMPERATURE MEASUREMENT	9
Calibration	9
Measurement	9
DATA COMMUNICATION	10
Recorder output	10
GLP-protocol	10
Digital output	11
Printer connection	11
Computer connection	11
TROUBLE SHOOTING	12
Simple test	12
Calibrate	12
Reset	12

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This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

Introduction

The normal operating temperature should be between 4° and 40°C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

If it is equipped with rechargeable batteries, the instrument should be connected permanently to the mains in order to keep the batteries in good condition. Do not store the instrument for longer than 1 month without recharging the batteries. The batteries can be replaced by opening the bottom compartment of the cabinet.

CONSORT nv
Parklaan 36
B2300 Turnhout
Belgium

Tel (++32)(14)41 12 79
Fax (++32)(14)42 91 79

Manufacturer

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

Warranty

CONSORT is not liable for consequential damages arising out of the use or handling of its products.

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply and **NOT** to send the following items unless they are suspect:

Servicing

Manual
Cables
Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

Ranges**Conductivity**

cc = 0.1 cm⁻¹ : 0.001 μS/cm ... 20 mS/cm
cc = 1 cm⁻¹ : 0.01 μS/cm ... 200 mS/cm
cc = 10 cm⁻¹ : 0.1 μS/cm ... 2000 mS/cm

Resistivity

cc = 0.1 cm⁻¹ : 100 Ω.cm ... 200 MΩ.cm
cc = 1 cm⁻¹ : 10 Ω.cm ... 20 MΩ.cm
cc = 10 cm⁻¹ : 1 Ω.cm ... 2 MΩ.cm

Salinity

cc = 0.1 cm⁻¹ : 0.001 ppm ... 2 %
cc = 1 cm⁻¹ : 0.01 ppm ... 10 %
cc = 10 cm⁻¹ : 0.1 ppm ... 10 %

°C 0...100°C

Specifications**Resolution**

0.1°C, 0.001 μS/cm, 1 Ω.cm, 0.01 ppm

Inputs

BNC input for conductivity cell
BANANA input for a Pt1000 temperature probe

Temp. Comp.

automatic with Pt1000 or manual

Digital output

programmable RS232, 2400 b/s, for bi-directional communication with a computer or printer

Analogue output

0...2 V, ca 8 kΩ, 8 bit D/A

Display

8 digit/character, 12 mm LCD display, shows readings, °C & messages

Keys

10 tactile membrane keys

Ambient temp.

4...40°C

Rel. Humidity

0...90 % (non-condensing!)

Power supply

210-250 V~ (* 105-125 V~), 50/60 Hz, max. 2 VA, & 4 NiCd batteries, size AA

Cabinet

IP65 cabinet

Dimensions

252 x 121 x 50 mm

Weight

600 g

Keyboard

- MODE** = Selects all modes or escapes from error traps, calibration procedures, etc..by returning to the original mode.
- CAL** = Starts or proceeds a calibration or a function.
- ▲/▼** = Button for entering a value or for selecting a function.
- °C** = Reads temperature, when a Pt1000 is present, or manual temperature compensation.
- HOLD** = Holds display when measuring.
- RES** = Changes resolution/units of the display.
- PRINT** = Displayed value is printed through the RS232 output.
- ON/OFF** = Switches the instrument on or off.

Each time the instrument is switched on, it will check briefly its internal calibrations while showing all functions of the display and the software-version. An error message appears when something is wrong.

Error codes

- [Or. EC]** = Overrange Conductivity (e.g. short circuited electrode cable).
- [Or. OhM]** = Overrange Resistivity (e.g. interrupted electrode cable).
- [Or. tdS]** = Overrange Salinity (e.g. short circuited electrode cable).
- [Or. °C]** = Overrange °C (measured degrees out of range).
- [no dAtA]** = Data-logging memory empty.
- [not.CAL]** = Calibration procedure interrupted (the previous calibration remains valid!).
- [not drY]** = Cell too wet for automatic capacitive compensation.
- [diLutE]** = Solution too concentrated for a TC plotting.
- [Err.cc]** = Cell constant not within 70...130% of the nominal value.
- [Err.CAL]** = Calibration error (check electrodes and standard solutions).
- [Err.Stb]** = Electrode not stable.
- [Err.EC]** = No cell connected during a TC plotting.
- [Err.Prt]** = Communication error with the attached printer.
- [Err.bAt]** = Battery charger defective (disconnect the AC-adapter immediately from the mains, remove the batteries and have your instrument serviced!).
- [Err.COM]** = General communication error.
- [Err.MEM]** = Fatal memory error (all factory settings are cleared. Have your instrument serviced!).
- [ILLEGAL]** = An unauthorised person has failed to repair the unit.
Warranty is elapsed from now on!

AC Adapter

Remove the protective silicone stopper from the REF socket and connect the jack of a suitable AC adapter (model **A4070** for 230 V~ or model **A4071** for 115 V~) to the socket for recharging the NiCd batteries. **Avoid to use an AC adapter without NiCd batteries being inserted in the instrument!** For field work, remove the adapter and reinsert the protective silicone stopper into the socket.

SET-UP

The built-in beeper can be programmed to sound each time a key-path is pressed.

Beeper

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with \blacktriangle or \blacktriangledown between [bEEP on] or [bEEP OFF]. Decide whether or not the beeper should sound each time the keyboard is touched.
3. Press **CAL** to continue or press **MODE** to cancel.

An audible alert can be given when the displayed values stray outside a programmable high- and low-level.

Alarm

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with \blacktriangle or \blacktriangledown between [ALrM OFF], [ALrM EC], or [ALrM tdS]. Press **CAL**. Select [ALrM OFF] if no alarm is required.
3. The display shows e.g. [L=23.0MS] while the units are blinking. Select the desired range with \blacktriangle or \blacktriangledown and press **CAL**.
4. The display shows e.g. [L=23.0MS] while [=] is blinking. Select the desired minimum level with \blacktriangle or \blacktriangledown and press **CAL**. Each time the measurements are below this level, the beeper will sound an alarm.
5. The display shows e.g. [H=57MS] while the units are blinking. Select the desired range with \blacktriangle or \blacktriangledown and press **CAL**.
6. The display shows e.g. [H=57MS] while [=] is blinking. Select the desired maximum level with \blacktriangle or \blacktriangledown and press **CAL**. Each time the measurements are above this level, the beeper will sound an alarm.
7. Press **CAL** to continue or press **MODE** to cancel.

The measured values can be printed automatically at timed intervals or manually on operator command.

Printer

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Pt=120]. Select the desired time interval (0...9999 s) between the transmitted data to a printer, with \blacktriangle or \blacktriangledown . Pre-set to zero if no automatic printing is required.
3. Press **CAL** to continue or press **MODE** to cancel.

The instrument can be identified by e.g. a computer when a specific number is allocated to it.

Identification number

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Id.no=5]. Select the desired identification number (0...999) for your meter with \blacktriangle or \blacktriangledown . Pre-set to zero if this number should not be mentioned while printing.
3. Press **CAL** to continue or press **MODE** to cancel.

A private code can be programmed to avoid undesired access to the instrument.

Secret code

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows [Code on], [Code OFF], or [Code CAL]. Decide, with \blacktriangle or \blacktriangledown , whether the access to all future calibrations should be by entering a code (on) or not (off). Press **CAL** to continue or press **MODE** to cancel. Select [Code CAL] to enter your personal code.
3. The display shows [E=-----] while [=] is blinking. Enter your secret sequence of 5 keys to proceed. Press **CAL** to continue or press **MODE** to cancel.

Permits to insatall the temperature compensation and the reference temperature.

Temperature compensation

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows [AtC=on], [MtC=on] or [AtC=OFF], [MtC=OFF]. Switch the temperature compensation on or off (for absolute measurements) with \blacktriangle or \blacktriangledown and press **CAL** to continue or **MODE** to return.
3. The display shows [rEF=25°C] or [rEF=20°C]. Select the temperature to which all future conductivity measurements will be referred with \blacktriangle or \blacktriangledown and press **CAL** to continue or **MODE** to return.

DATA-ACQUISITION

Up to 100 values can be stored in the non-volatile memory of the instrument.

- [no LOG] : No data-logging.
[SEt LOG] : Permits to pre-set the data-logging parameters.
[LSt LOG] : Permits to view the stored values on the display.
[Prt LOG] : Permits to send the stored values to a printer or computer.
[Cnt LOG] : Permits to continue with a next series of measurements, without erasing the previous data-logging.

Set log: Procedure to pre-set the data-logging parameters.

Logging data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [SEt LOG] with \blacktriangle or \blacktriangledown . Select [no LOG] if no data-logging is required and press **MODE** to return to the original mode.
3. Press **CAL** until you can choose with \blacktriangle or \blacktriangledown between [LOG EC], [LOG OhM] or [LOG tdS]. Select the desired mode and press **CAL**.
4. The display shows e.g. [dt=100]. Select the desired time interval (0...9999 s) between the data-logging, with \blacktriangle or \blacktriangledown and press **CAL**. Pre-set to zero for manual data-logging, each time **PRINT** is pressed (the next points 5 to 7 are skipped and press **CAL**).
5. The display shows e.g. [n=40]. Select the desired number of values (1...100) to be data-logged with \blacktriangle or \blacktriangledown and press **CAL**.

6. The display shows e.g. [d=31.12.91] while the year is blinking. Program the desired starting date with \blacktriangle or \blacktriangledown and press **CAL**. Program similarly the month and the day.
7. The display shows e.g. [t=23.59] while the minutes are blinking. Program the desired starting time with \blacktriangle or \blacktriangledown and press **CAL**. Program similarly the hours.
8. The display shows [StArt ?]. Press **CAL** to start the data-logging according the previous settings (a blinking [L] appears) or press **MODE** to cancel. After the last value has been catalogued the display will show [End LOG]. When manual data-logging has been selected, press **PRINT** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [LOG 7]. Press **MODE** to cancel.

[Cnt LOG]: Permits to continue with a next series of measurements, without erasing the previous data-logging

Adding data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Cnt LOG] with \blacktriangle or \blacktriangledown to continue with a next series of measurements, without erasing the previous data-logging, provided the memory has enough capacity left and press **CAL**.

[Lst LOG]: permits to view the stored values on the display.

Recalling data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Lst LOG] with \blacktriangle or \blacktriangledown and press **CAL**
3. The display shows the first data-logged value in memory (a blinking [n] appears). Press \blacktriangle or \blacktriangledown to view all other values.
4. While pressing **HOLD** the corresponding date is shown (not relevant when a manual or continued data-logging has been performed).
5. While pressing **RES** the corresponding time shown (or: series + log-number, e.g. [2-017], when a continued data-logging has been performed).
6. While pressing **°C** the corresponding temperature is shown.
7. Press **PRINT** to send the recalled data to a printer.
8. Press **MODE** to return to cancel.

[Prt LOG]: Permits to send the stored values to a printer or computer.

Printing data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [Prt LOG] with \blacktriangle or \blacktriangledown and press **CAL**
3. The display shows [PrintinG] while all data-logged values, including corresponding dates and times, are sent to the RS232 output in the following format, compatible with EXCEL™, LOTUS 123™,... The columns are separated by a TAB character.

Log#	VALUE	UNIT	C	H	D
001	16.49E-03	S/cm	23.7	13:19:57	29/01/95
002	15.05E-03	S/cm	23.7	13:20:57	29/01/95
003	13.39E-03	S/cm	23.8	13:21:57	29/01/95

The measuring cell should be connected to the coaxial INPUT connector. Automatic temperature compensation and temperature measurements are possible by plugging a Pt1000 thermocompensator into the °C terminals. You can also use a combination conductivity electrode with built in Pt1000. Its banana plug should be inserted in the upper °C terminal. Without compensator, the manual temperature compensation is automatically switched on.

CONDUCTIVITY MEASUREMENT

A conductivity cell is active and stable only after wetting! For this purpose it must be immersed for **at least one hour** in distilled water. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

A polluted cell may be cleaned with a soft detergent or diluted nitric acid. Greasy substances may be removed with acetone or alcohol (never do this with plastic electrodes!).

Maintenance

The cell constant depends on the physical dimensions of the platinum plates. They are coated with a spongy black platinum deposit. This increases greatly the effective surface and reduces polarising effects. Since the coating of platinum black is quite resistant to chemical action, the cell constant remains fairly constant during the electrode life time.

Calibration

1. Select the conductivity range by pressing **MODE** (the display shows [μS] or [MS]) and press **CAL**.
2. The display will show the cell constant (e.g. [cc=1.073]). Press **CAL** to start the automatic determination procedure.
3. The display shows one of the KCl standard solutions in memory [St=0.01M] (1413 μS/cm), [St=0.1M] (12.88 mS/cm), [St=1M] (111.8 mS/cm) or the manual input [St=MAN]. Select the proper standard with ▲ or ▼.
4. This step is **skipped** unless the manual input has been chosen. The display shows e.g. [MS=10.00] while the units are blinking. Choose with ▲ or ▼ between 1.000 (units), 10.00 (tens) or 100.0 (hundreds) the desired calibration range and press **CAL** again. The display shows e.g. [MS=1.234] while [=] is blinking. Select the desired standard value with ▲ or ▼.
5. After rinsing the cell several times with the standard solution, immerse it the standard solution. The solution temperature is not so critical but it should lie between 0°C and 30°C. If the standard has a different temperature press °C here (not necessary when using a Pt1000). Press **CAL** to continue.
6. The display will show the computed cell constant e.g. [cc1=1.045] and will calibrate automatically when readings are stable ([=] stops blinking).

Temperature compensation

Each solution has its own temperature coefficient, which also varies with the temperature. Therefore it is impossible to achieve an exact temperature compensation over a wide span of temperatures with a conventional conductometer. However, this model features a unique microprocessor controlled system, which helps you to completely overcome the problem of non linear temperature compensation curves for most of the species to be measured. There is no need to keep sample temperatures close to each other, since you can plot special temperature curves from up to 10 different solutions.

1. Press **MODE** until the display will show e.g. [tc nACI], [tc07 PrG] or [tc07 CLr] depending on temperature curve 07 being programmed previously or being cleared. Remember that [tc nACI] is the fixed NaCl temperature curve which can never be altered. Select the desired temperature curve with ▲ or ▼ and press **MODE** to proceed or **CAL** to start the automatic tracing.
2. Prepare the sample solution to be measured in a thermostatic bath with magnetic stirrer. Its temperature should be about 5°C lower than the lowest value you may possibly need during future measurements, but at least lower than 10°C. The display shows e.g. [tc07] [8.9] which means temperature curve 07 at 8.9°C. After rinsing the electrodes with the solution to be measured, immerse them in that solution. Wait a while to enable the electrodes to adapt themselves to the low temperature. Press **CAL** again to start the automatic plotting procedure.
3. Slowly increase the temperature (max. 0.1°C/s) of the thermostatic bath. You can monitor the temperature on the display while a blinking [c], indicating that plotting has started, appears. When the temperature is about 5°C higher than the highest value you need for future measurements (min. 40°C required), press **CAL** to stop the temperature tracing. The display shows shortly [WAlt] while the curve is being computed. After the display shows e.g. [tc07 PrG], select any mode to proceed.

- * After a temperature curve has been selected and **PRINT** is pressed, the instrument transmits the whole curve to the digital output. In the meantime the display shows shortly [PrintinG].
- * If you have traced the temperature compensation in a limited range (e.g. from 10° to 40°C), the curve is supposed to go on linearly outside these limits.

1. Verify if the attached cell is completely dry.
2. Select the conductivity range by pressing **MODE** (the display shows [μS] or [MS]) and press **CAL**.
3. The display will show the cell constant (e.g. [cc=1.073]), [COMP.on] or [COMP.OFF]. Select [COMP.on] with ▲ or ▼ and press **CAL** to start the automatic capacitance compensation. The display will show [WAlt] for a while.

Capacitance compensation

Conductivity measurement

1. Select the proper temperature compensation with **MODE**. If you ignore this step, the instrument will automatically assume the last selected temperature curve
2. Select the conductivity range by pressing **MODE** (the display shows [μ S] or [MS]).
3. Rinse the electrodes with the sample and immerse them in it.
4. Read the display (remember: compensated to the reference temperature!). The autoranging feature gives you the largest possible number of digits + their unit.
5. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

- * Disconnect the AC-adapter to increase accuracy in the lower conductivity range!
- * See page 49 for the standards in memory.
- * A blinking decimal point warns you for unstable measurements. Wait to read the display!
- * Press **RES** to switch between ppm (salinity), S/cm (conductivity) or Ω .cm (resistivity).
- * When starting a titration, press ∇ to lock the actual range and avoid cross-over errors due to differences in the measuring frequency. Press **MODE** to return to the normal measuring mode.
- * Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.

Salinity measurement

1. Select the conductivity range by pressing **MODE** (the display shows [μ S] or [MS]) and press **RES** until the display shows ppm (salinity range).
2. Rinse the electrodes with the sample and immerse them in it.
3. Read the display. The autoranging feature gives you the largest possible number of digits + their unit.
4. Rinse the cell always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).

Relative measurements

1. While in the conductivity or salinity mode, press \wedge . At that moment the instrument stores the actual value and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
2. To cancel relative readings, press **MODE** and you will be in the normal measuring mode again.

TEMPERATURE MEASUREMENT

1. Press **CAL** to calibrate the $^{\circ}$ C scale with any reference. Immerse the Pt1000 in a solution of known temperature.
2. The display shows e.g. [$^{\circ}$ C=22.3]. Calibrate to the proper value with \wedge or ∇ and press **CAL** again or press **RES** to reset the absolute readings.

Calibration

1. Read temperature by pressing °C.
2. Without Pt1000, adjust the manual temperature compensation with ▲ or ▼ and proceed by pressing **MODE**.

Measurement

DATA COMMUNICATION

A recorder can be connected to the red (+) and the black (-) terminals. The 0 to 2 V output range corresponds with 0 to 2000 digits. Use only laboratory recorders with a high input impedance!

Recorder output

1. Select the conductivity range by pressing **MODE**.
2. Press **CAL** to display the cell constant, e.g. [cc=1.073].
3. Press **PRINT** to print a complete calibration report, e.g.:
4. Press **MODE** to return to the normal measuring mode.

GLP-protocol

```

CONSORT Conductometer K612
-----

SETTINGS
-----
Identification No      : 004
Software Version      : 3.0
User Code              : on
Titration Lock        : off
Rel. Conductivity     : off
Cap. Compensation     : off

ATC      (Pt1000)     : on
Calibr. Temp. (C)    : 23.4
Ref. Temp. (C)      : 25.0

Alarm EC              : on
Alarm High (S/cm)    : 85.7E-03
Alarm Low (S/cm)     : 30.5E-03

Temp. Comp.          : NaCl

Batt.Capacity (%)    : 87

CALIBRATION
-----
Ref. Temp. (C)      : 25
Standard (S/cm)     : 1413E-06
Cell Const (cm-1)   : 1.073
Response Time (s)   : 13

STATISTICS
-----
Cell Const (cm-1)   : -0.026
Average Resp. (s)   : +1

```

Digital output

A standard RS232 output terminal (DP9) is provided for interfacing the instrument with a printer or computer. Data is sent in the ASCII code at a BAUD rate of 2400 b/s (8 bit, no parity, 1 start and 2 stopbits, automatic line feed ON, data transfer protocol Xon/Xoff).

Serial port pinout specifications:

pin 1 : RLSD, received line signal detector
pin 2 : TxD, transmit data
pin 3 : RxD, receive data
pin 4 : not connected
pin 5 : Gnd, signal ground
pin 6 : DSR, data set ready
pin 7 : not connected
pin 8 : CTS, clear to send
pin 9 : not connected

Interfacing with a printer:

Press **PRINT** each time you want to print the measured values. The columns are separated by a TAB character. Run the set-up procedure (page for automatic printing at timed intervals, 'Printer' page 4, while you are measuring. The **PRINT** keypath does not function in this case. It is also disabled during an automatic data-logging., e.g.:

Log#	VALUE	UNIT	C	H	D
001	16.49E-03	S/cm	23.7	13:19:57	29/01/95
002	15.05E-03	S/cm	23.7	13:20:57	29/01/95
003	13.39E-03	S/cm	23.8	13:21:57	29/01/95

Printer connection

Using WINDOWS™ to collect data:

1. Connect the computer with the RS232 terminal, using the optional AK2310 cable (eventually with adapter AB3190 for connection with a 25 pole RS232 terminal).
2. Start **WINDOWS™** and select *Terminal* in the *Accessories Group*.
3. Choose *Receive Text File* in the *Transfer Menu* and create a file name (TXT format) for the data to receive.
4. Select the *Settings Menu* and make sure the following parameters are programmed:

Baud Rate : 2400
Data bits : 8
Stop bits : 2
Parity : None
Flow control : Xon/Xoff

5. Start the conductometer in the desired measuring mode for manual or automatic data collection (see 'Printer' page 4 for more details).
6. When finished, open the created file with a standard spreadsheet (e.g. **EXCEL™**) or wordprocessor (e.g. **WORD™**) for further processing.

Computer connection

Interfacing with a computer:

Send a single character to the instrument and it will execute the corresponding command as follows:

- 1 = press **MODE**.
- 2 = press **CAL**.
- 3 = press **▲**.
- 4 = press **▼**.
- 5 = press **SET**.
- 6 = press **HOLD**.
- 7 = press **RES**.
- 8 = press **PRINT**.
- ? = send display to computer.
- + = keyboard on (manual operation remains possible).
- = keyboard off (manual operation disabled).

Follow exactly the same measuring or calibration procedures, as described in this manual, to include any desired command in your computer data-acquisition program.

This example continuously prints the measured values (pH and temperature) on the computer screen:

10	ON ERROR GOTO 70	'makes sure no error occurs
20	OPEN "COM1:2400,N,8,2" AS #1	'prepares computer input
30	PRINT #1,"8"	' PRINT button pressed
40	INPUT #1,V\$	'measured values transmitted
50	PRINT V\$	'received values printed
60	GOTO 30	'repeat procedure
70	RUN	

TROUBLE SHOOTING

Try this procedure if you suspect the electronic performance of the instrument.

1. Leave the conductivity input open.
2. Switch the instrument ON.
3. Select the conductivity range by pressing **MODE**. The display should show $[0.00\mu\text{S}] \pm 0.01 \mu\text{S/cm}$. Run the **Calibrate** procedure if erroneous values are read.

This procedure permits to calibrate automatically all amplifiers of the instrument.

1. Leave the conductivity input open.
2. Switch the instrument ON while holding **CAL** pressed.
3. The display shows [SELFCAL] for a while and returns to the measurements.

This procedure resets the instrument to the original factory settings. All personal memorised data (including access code) will be erased!

1. Switch the instrument ON while holding **MODE** pressed.
2. The display shows [rESET] for a while and returns to the measurements.

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