
Consort

**MANUAL
HANDLEIDING
MODE D'EMPLOI
ANLEITUNG**

**T8710
T8720**

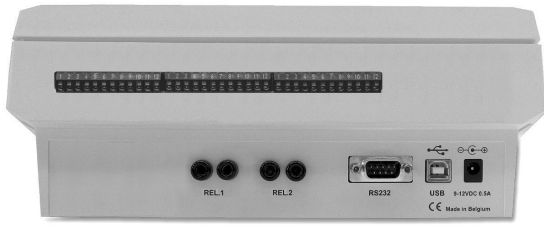April 2013

T8710

16 channels for thermocouples type J, K, T or E

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General informations

Introduction

This instrument is manufactured with the latest technology and needs no particular maintenance. **Consort** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 0° and 40° C. Never use the instrument in a room with high humidity (>95 %) or at very low temperatures (condensation water!).

Manufacturer

Consort bvba
Parklaan 36
B2300 Turnhout
Belgium

Tel (+32)(14)41 12 79
Fax (+32)(14)42 91 79
E-mail: info@consort.be

Warranty

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **Consort** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **Consort** before returning any product for warranty repair on a freight prepaid basis!

Consort is not liable for consequential damages arising out of the use or handling of its products.

Servicing

In the event of this instrument being returned for servicing, the owner is requested **NOT** to send the following items unless they are suspect:

Manual

Cables

Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **Consort** dealer.

AC adaptor

Connect the mains adapter to the DC socket. Do not hold the adapter by wet hand.

Keyboard

MODE = Selects the settings or escapes from error traps, calibration procedures, etc..by returning to the original mode.

↔↑↓ = Button for entering a value, for selecting a function or a channel.

CAL = Starts or proceeds a calibration or a function.

STORE = Stores the displayed value or send it to a computer.

HOLD = Holds display when measuring.

ON/OFF = Switches the instrument on or off.

Inputs

T8710

Connect the temperature probes with the terminal blocks and insert them in the instrument. The channel numbers start from the left to the right (backview). The even numbers are the negative poles. The uneven numbers are the positive poles.

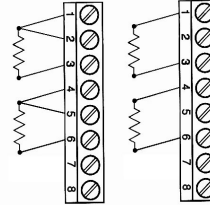
T8720

Connect the temperature probes with the terminal blocks and insert them in the instrument. The channel numbers start from the left to the right (backview).

2-wire Pt100: connect to clamps 1 and 3, etc.

3-wire Pt100: connect to clamps 1+2 and 3, etc.

4-wire Pt100: connect to clamps 1+2 and 3+3, etc.

Outputs

A standard USB or Ethernet port is provided for interfacing the instrument with a computer or network.

A standard RS232 port is provided for interfacing the instrument with a printer. Data is sent in the ASCII code at a BAUD rate of 300...19200 bps (8 bit, no parity, 1 stopbit).

- The RS232 port is automatically switched off when a USB cable is connected.

Language

1. Press **MODE**.
2. Select [SYSTEM] and press **CAL**.
3. Select [LANGUAGE] and press **CAL**.
4. Select the desired language and press **CAL**.

Contrast

1. Press **MODE**.
2. Select [SYSTEM] and press **CAL**.
3. Select [CONTRAST] and press **CAL**.
4. Adjust the contrast and press **CAL**.

Date and time

1. Press **MODE**.
2. Select [SYSTEM] and press **CAL**.
3. Select [DATE/TIME] and press **CAL**.
4. Select date or time and press **CAL** to adjust.

Password

1. Press **MODE**.
2. Select [PASSWORD] and press **CAL**.
3. A private code can be programmed to avoid undesired access to the instrument. Select [YES] and press **CAL** to enter your secret sequence of 5 keys.

Channels

1. Press **MODE**.
2. Select [CHANNELS] and press **CAL**.
3. Select [ACTIVATE] and press **CAL**.
4. Select the desired channel and press **CAL** to activate or deactivate.
5. Select [THERMOCOUPLE] and press **CAL**. (*T8710 only*)
6. Select the desired channel and press **CAL**. (*T8710 only*)
7. Select the type of thermocouple and press **CAL**. (*T8710 only*)
8. Select [UNIT] and press **CAL**.
9. Select the desired unit and press **CAL**.
10. Select [RESOLUTION] and press **CAL**.
11. Select the desired measurement resolution and press **CAL**.

Stand-by

1. Press **MODE**.
2. Select [SYSTEM] and press **CAL**.
3. Select [STAND-BY] and press **CAL**.
4. Select [YES] and press **CAL** to restart automatically after a power failure.

Relays

1. Press **MODE**.
2. Select [RELAYS] and press **CAL**.
3. Select [ACTION] for relay-1 and press **CAL**.
4. Select [OPEN] [CLOSE] or [NONE] and press **CAL**.
5. Select [WHEN] for relay-1 and press **CAL**.
6. Select when the relay should be closed and press **CAL**.:
 - [ALARM] maximum or minimum level is exceeded.
 - [ALARM>] maximum level is exceeded.
 - [ALARM<] minimum level is exceeded.
 - [STAND-BY] instrument is in stand-by.
 - [ACTIVE] instrument is measuring.
 - [INACTIVE] instrument is not measuring.
7. Proceed for relay-2 accordingly.
 - *Specifications of the relays: max. 50 V, 200 mA, AC/DC*

Alarm

1. Select the menu by pressing **MODE**.
2. Select [ALARM] and press **CAL**.
3. Select [BEEPER] and press **CAL** to activate or deactivate.
4. Select the desired channel and press **CAL**.
5. Select [ACTIVATE] and press **CAL** to activate or deactivate.
6. Select [TEMP.>] and press **CAL**.
7. Select the desired maximum level and press **CAL**. Each time the measurements are above this level, the beeper will sound an alarm.
8. Select [TEMP.<] and press **CAL**.
9. Select the desired minimum level and press **CAL**. Each time the measurements are below this level, the beeper will sound an alarm.
 - *Alarm hysteresis = 1 °C*

Calibration

1. Verify that the instrument is switched on in a stable temperature environment.
2. Press **CAL** while measuring.
3. Select [CHANNEL] and press **CAL**.
4. Select the desired channel and press **CAL**.
5. Select [METHOD] and press **CAL**.
6. Select the desired calibration method and press **CAL**.
7. Select [TYPE] and press **CAL**.
8. Select the 1- or 2-point calibration and press **CAL**.
9. Select [CALIBRATE] and press **CAL**. Follow the instructions on the screen.

Measurement

1. Verify that the instrument is switched on in a stable temperature environment. The display immediately shows the measured value according to the last calibration.
 - Press **↔** to change between the channels.
 - Press **↑** to display the maximum temperature ever measured by the selected channel (after switching on). Press **↑** again to return to the normal measurements. Press **CAL** to reset the minimum/maximum memory.
 - Press **↓** to display the minimum temperature ever measured by the selected channel (after switching on). Press **↓** again to return to the normal measurements. Press **CAL** to reset the minimum/maximum memory.
 - Any alarm will sound an alert and close a relay when programmed for alarm. Press **MODE** to stop the beeper.

Alternating measurements

1. Select the menu by pressing **MODE**.
2. Select [ALTERNATE] and press **CAL** to activate or deactivate. The display alternates between the selected modes with an interval of ca. 5 s.
3. Press **MODE** to return to the measurements.

Good Laboratory Practice

1. Press **CAL** while measuring.
2. Select [GLP] and press **CAL**.
3. Browse with **←→↑↓** to show a complete calibration report. Press **STORE** to print the report.

Data Storage in the internal memory

1. Select the menu by pressing **MODE**.
2. Select [DATALOGGER] and press **CAL**.

Start the data-logging

1. Select [STORE] and press **CAL**.
2. Select [ACTIVATE] and press **CAL**.
3. Select [YES] or [NO] and press **CAL**.
4. Select [INTERVAL] and press **CAL**.
5. Select the desired time interval between the data-logging (or MANUAL) and press **CAL**.
6. Select [ALARM] and press **CAL**.
7. Select the desired time interval between the data-logging (or MANUAL), when the alarm is activated, and press **CAL**.
8. Select [NUMBER] and press **CAL**.
9. Select the desired number of values to be data-logged and press **CAL**.
10. Press **MODE** to return to the measurements and to start the data-logging according to the previous settings. When manual data-logging has been selected, press **STORE** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [#00027].

View the stored values on the display

1. Select [RECALL] and press **CAL**.
2. Select [TABLE] or [PRINT] to display or print the stored data and press **CAL** to continue. Follow the instructions on the screen.
3. In case [TABLE] have been selected press **CAL** to show the options.
4. Select [LOG#] and press **CAL**. Select the number of the first line to be shown on the display and press **CAL**.
5. Select [SHOW] and press **CAL**. Select [Log#] (line numbers) or [Date] to be shown on the display and press **CAL**.

Erase the stored values

1. Select [ERASE] and press **CAL**. Follow the instructions on the screen.

External storage in a computer using USB or RS232

1. Select the menu by pressing **MODE**.
 2. Select [USB/RS232] and press **CAL**.
 3. Select [BAUDRATE] and press **CAL**.
 4. Select the desired baudrate and press **CAL**.
 5. Select [INTERVAL] and press **CAL**.
 6. Select the desired interval between the transmitted data and press **CAL**. Pre-set to zero if no automatic transmitting is required.
 7. Select [ALARM] and press **CAL**.
 8. Select the desired interval between the transmitted data, when the alarm is activated, and press **CAL**. Pre-set to zero if no automatic transmitting is required.
 9. Select [IDENTIF. No.] and press **CAL**.
 10. Enter an identification number for the transmitted data and press **CAL**.
- You can download free software and a suitable USB driver from www.consort.be.

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Specifications		T8720
TEMPERATURE	Range, Pt100	-200...850°C (-328...1562°F)
	Resolution	0.1°C (0.1°F)
	Accuracy	0.1% ±0.3°C (±0.5°F)
	Calibration	1...2 points
	Channels	12
INPUTS		terminal blocks for 12 Pt100 probes
CALIBRATION	GLP Protocol	✓
DISPLAY	LCD	128x64 pixels
	Alternating	✓
	Hold Function	✓
	Selectable Resolution	✓
	Real Time Clock	✓
	Languages	EN, NL, FR, DE
COMMUNICATION	Interface with computer	USB
	RS232, baud rate	300...115200 b/s
	Printer	✓
DATA-LOGGING	Values	18000 + date/time
	Manual	✓
	Timed	✓
	Interval	4...9999 s
ALARM	Relays	2, max. 50 V/200 mA
SECURITY	Identification Number	✓
	Password Protection	✓
AMBIENT CONDITIONS	Temperature	0...40°C
	Humidity	0...95%, non condensing
POWER SUPPLY	Mains	90...250 VAC, 50/60 Hz
DIMENSIONS	WxDxH	26x18x9 cm
WEIGHT	Meter	1 kg

DECLARATION OF CONFORMITY

We declare under our sole responsibility that the product

**Multi-channel analysers
content of the type numbers
T8710, T8720**

to which this declaration relates is in conformity
with the following standards

LOW VOLTAGE DIRECTIVE 2006/95/EC

EN61010-1

EMC DIRECTIVE 2004/108/EC

EN61326-1

Consort bvba

Parklaan 36, B-2300 Turnhout, Belgium

Tel: (+32) (0)14 41 12 79

Fax: (+32) (0)14 42 91 79

Sales: sales@consort.be

The electro-magnetic susceptibility has been chosen at a level that gains proper operation in residential areas, on business and light industrial premises and on small-scale enterprises, inside as well as outside of buildings. All places of operation are characterised by their connection to the public low voltage power supply system.

WARRANTY CERTIFICATE

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory.

Accessories and breakable items such as electrodes are not warranted unless proven to be defective before shipment.

The original purchase order numbers, Consort invoice numbers and serial numbers of the products must be provided.

CONSORT will repair all defective equipment returned to it during the warranty period without charge (CIF Turnhout prepaid by sender), provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures described in the instruction manual and when not having been subject to accident, alteration, misuse or abuse.

If the products have been used with or have come into contact with fluids, an MSDS (material safety data sheet) must be supplied prior to issuing a return authorisation.

A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

Consort bvba

Parklaan 36 • B-2300 Turnhout • Belgium

Tel: (+32) (0)14 41 12 79

Fax: (+32) (0)14 42 91 79

Sales: sales@consort.be

Support: support@consort.be

Information: info@consort.be