# CONSORT

---

MANUAL

HANDLEIDING

MODE D'EMPLOI

ANLEITUNG

**Z621**

---

January 1997

## **Table of contents**

Introduction .....	1
Manufacturer .....	1
Warranty .....	1
Servicing .....	1
Specifications .....	2
Keyboard .....	2
Error codes .....	3
<b>SET-UP .....</b>	<b>3</b>
Beeper .....	3
Alarm .....	3
Printer .....	4
Identification number .....	4
Secret code .....	4
<b>DATA-ACQUISITION .....</b>	<b>4</b>
Logging data .....	5
Adding data .....	5
Recalling data .....	5
Printing data .....	6
Inputs .....	6
<b>MEASUREMENTS .....</b>	<b>6</b>
Theory .....	6
Maintenance .....	7
Measurements	
with 1-point calibration .....	7
with 2-point calibration .....	8
Relative measurement .....	9
Temperature measurement .....	9
<b>DATA COMMUNICATION .....</b>	<b>9</b>
Recorder output .....	9
GLP-protocol .....	9
Digital output .....	10
Printer connection .....	11
Computer connection .....	11
<b>TROUBLE SHOOTING .....</b>	<b>12</b>
Calibrate .....	12
Reset .....	12

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This instrument is manufactured with the latest technology and needs no particular maintenance. **CONSORT** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

## ***Introduction***

The normal operating temperature should be between 4° and 40°C. Never store the instrument in a room with high humidity or at very low temperatures (condensation water!).

If it is equipped with rechargeable batteries, the instrument should be connected permanently to the mains in order to keep the batteries in good condition. Do not store the instrument for longer than 1 month without recharging the batteries. The batteries can be replaced by opening the bottom compartment of the cabinet.

**CONSORT nv**  
Parklaan 36  
B2300 Turnhout  
Belgium

Tel (++32)(14)41 12 79  
Fax (++32)(14)42 91 79

## ***Manufacturer***

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory. **CONSORT** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

## ***Warranty***

**CONSORT** is not liable for consequential damages arising out of the use or handling of its products.

In the event of this instrument being returned for servicing, the owner is requested to remove the power supply lead and **NOT** to send the following items unless they are suspect:

## ***Servicing***

*Manual*  
*Cables*  
*Accessories*

If serious malfunctioning occurs, stop using the unit immediately and consult your local **CONSORT** dealer.

## Specifications

<b>Ranges</b>	<b>O<sub>2</sub></b> 0 ... 60 ppm <b>Saturation</b> 0 ... 600 % <b>°C</b> 0 ... 50°C
<b>Accuracy</b>	0.01 ppm/ 0.1 %/ 0.1°C or 1% of reading
<b>Inputs</b>	BNC + BANANA input for oxygen electrode
<b>Temp. comp.</b>	automatic with NTC (100 kΩ), built-in O <sub>2</sub> -probe
<b>Digital output</b>	programmable RS232, 2400 b/s, for bi-directional communication with a computer or printer
<b>Analogue output</b>	0 ... 2 V, ca 8 kΩ, 8 bit D/A
<b>Display</b>	8 digit/character, 12 mm LCD display, shows readings, °C & messages
<b>Keys</b>	10 tactile membrane keys
<b>Ambient temp.</b>	4 ... 40°C
<b>Rel. Humidity</b>	0 ... 90 % (non-condensing!)
<b>Power supply</b>	210-250 V~ (* 105-125 V~), 50/60 Hz, max. 2 VA, & 4 NiCd batteries, size AA
<b>Cabinet</b>	IP65 cabinet
<b>Dimensions</b>	252 x 121 x 50 mm
<b>Weight</b>	600 g

Each time the instrument is switched on, it will check briefly its internal calibrations while showing all functions of the display and the software-version. An error message appears when something is wrong.

<b>MODE</b>	=	Selects all modes (except temperature) or escapes from error traps, calibration procedures..., by returning to the original mode.
<b>°C</b>	=	Reads temperature.
<b>▲/▼</b>	=	Button for entering a value or for selecting a function.
<b>CAL</b>	=	Starts or proceeds a calibration or function.
<b>HOLD</b>	=	Holds the display when measuring.
<b>RES</b>	=	Changes the resolution of the display.
<b>PRINT</b>	=	Displayed value is printed through the RS232 output.
<b>ON/OFF</b>	=	Switches the instrument on or off.

## Keyboard

## **Error codes**

- [Or.PPM] = Overrange ppm (e.g. defective electrode).  
[Or.%] = Overrange % (e.g. defective electrode).  
[Or.°C] = Overrange °C (measured degrees out of range).  
[no dAtA] = Data-logging memory empty.  
[not.CAL] = Calibration procedure interrupted (the previous calibration remains valid!).  
[Err.°C] = Temperature probe not connected.  
[Err.Prb] = Zero-point too large (e.g. defective membrane).  
[Err.Stb] = Electrode not stable during a calibration.  
[Err.CAL] = Calibration error (check electrode & standard solutions).  
[Err.COM] = General communication error.  
[Err.Prt] = Communication error with the attached printer.  
[Err.bAt] = Battery charger defective (disconnect the AC-cord immediately from the mains, remove the batteries and have your instrument serviced!).  
[Err.MEM] = Fatal memory error (all factory settings are cleared. Have your instrument serviced!).  
[ILLEGAL] = An unauthorised person has failed to repair the unit.  
**Warranty is elapsed from now on!**

## **SET-UP**

*The built-in beeper can be programmed to sound each time a key-path is pressed.*

### **Beeper**

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [bEEP on] or [bEEP OFF]. Decide whether or not the beeper should sound each time the keyboard is touched.
3. Press **CAL** to continue or press **MODE** to cancel.

*An audible alert can be given when the displayed values stray outside a programmable high- and low-level.*

### **Alarm**

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [ALrM OFF], [ALrM PPM] or [ALrM %]. Select [ALrM OFF] if no alarm is required. Press **CAL**.
3. The display shows e.g. [L=4.00PPM] or [L=60.0%]. Select the desired minimum level with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Each time the measurements are below this level, the beeper will sound an alarm.
4. The display shows e.g. [H=6.00PPM] or [H=80.0%]. Select the desired maximum level with  $\blacktriangle$  or  $\blacktriangledown$ . Each time the measurements are above this level, the beeper will sound an alarm.
5. Press **CAL** to continue or press **MODE** to cancel.

## Printer

*The measured values can be printed automatically at timed intervals or manually on operator command.*

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Pt=120]. Select the desired time interval (0...9999 s) between the transmitted data to a printer, with  $\blacktriangle$  or  $\blacktriangledown$ . Pre-set to zero if no automatic printing is required.
3. Press **CAL** to continue or press **MODE** to cancel.

*The instrument can be identified by e.g. a computer when a specific number is allocated to it.*

## Identification number

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows e.g. [Id.no=5]. Select the desired identification number (0...999) for your meter with  $\blacktriangle$  or  $\blacktriangledown$ . Pre-set to zero if this number should not be mentioned while printing.
3. Press **CAL** to continue or press **MODE** to cancel.

*A private code can be programmed to avoid undesired access to the instrument.*

## Secret code

1. Select the set-up procedure by pressing **MODE**.
2. Press **CAL** until the display shows [Code on], [Code OFF], or [Code CAL]. Decide, with  $\blacktriangle$  or  $\blacktriangledown$ , whether the access to all future calibrations should be by entering a code (on) or not (off). Press **CAL** to continue or press **MODE** to cancel. Select [Code CAL] to enter your personal code.
3. The display shows [E=-----] while [=] is blinking. Enter your secret sequence of 5 keys to proceed. Press **CAL** to continue or press **MODE** to cancel.

# DATA-ACQUISITION

*Up to 100 values can be stored in the non-volatile memory of the instrument.*

- [no LOG] : No data-logging.
- [Set LOG] : Permits to pre-set the data-logging parameters.
- [LSt LOG] : Permits to view the stored values on the display.
- [Prt LOG] : Permits to send the stored values to a printer or computer.
- [Cnt LOG] : Permits to continue with a next series of measurements, without erasing the previous data-logging.

## Set log: Procedure to pre-set the data-logging parameters.

## Logging data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [SEt LOG] with  $\blacktriangle$  or  $\blacktriangledown$ . Select [no LOG] if no data-logging is required and press **MODE** to return to the original mode.
3. Press **CAL** until you can choose with  $\blacktriangle$  or  $\blacktriangledown$  between [LOG PPM] or [LOG %]. Select the desired mode and press **CAL**.
4. The display shows e.g. [dt=100]. Select the desired time interval (0...9999 s) between the data-logging, with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Pre-set to zero for manual data-logging, each time **PRINT** is pressed (the next points 5 to 7 are skipped and press **CAL**).
5. The display shows e.g. [n=40]. Select the desired number of values (1...100) to be data-logged with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**.
6. The display shows e.g. [d=31.12.91] while the year is blinking. Program the desired starting date with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Program similarly the month and the day.
7. The display shows e.g. [t=23.59] while the minutes are blinking. Program the desired starting time with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**. Program similarly the hours.
8. The display shows [StArt ?]. Press **CAL** to start the data-logging according the previous settings (a blinking [L] appears) or press **MODE** to cancel. After the last value has been catalogued the display will show [End LOG]. When manual data-logging has been selected, press **PRINT** to put a next measurement into memory. In the meantime the display shows the logging-number e.g. [LOG 7]. Press **MODE** to cancel.

## [Cnt LOG]: Permits to continue with a next series of measurements, without erasing the previous data-logging

## Adding data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Cnt LOG] with  $\blacktriangle$  or  $\blacktriangledown$  to continue with a next series of measurements, without erasing the previous data-logging, provided the memory has enough capacity left and press **CAL**.

## [Lst LOG]: permits to view the stored values on the display.

## Recalling data

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Select [Lst LOG] with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL**
3. The display shows the first data-logged value in memory (a blinking [n] appears). Press  $\blacktriangle$  or  $\blacktriangledown$  to view all other values.
4. While pressing **HOLD** the corresponding date is shown (not relevant when a manual or continued data-logging has been performed).
5. While pressing **RES** the corresponding time shown (or: series + log-number, e.g. [2-017], when a continued data-logging has been performed).
6. While pressing  $^{\circ}\text{C}$  the corresponding temperature is shown.
7. Press **PRINT** to send the recalled data to a printer.
8. Press **MODE** to return to cancel.

**[Prt LOG]:** Permits to send the stored values to a printer or computer.

## *Printing data*

1. Select the set-up procedure by pressing **MODE** and press **CAL**.
2. Choose [Prt LOG] with **▲** or **▼** and press **CAL**
3. The display shows [PrintinG] while all data-logged values, including corresponding dates and times, are sent to the RS232 output in the following format, compatible with EXCEL™, LOTUS 123™,... The columns are separated by a TAB character.

Log#	VALUE	UNIT	C	H	D
001	7.34	ppm	23.7	13:19:57	29/01/95
002	7.21	ppm	23.7	13:20:57	29/01/95
003	7.08	ppm	23.8	13:21:57	29/01/95
004	6.71	ppm	23.8	13:22:57	29/01/95
005	6.49	ppm	23.9	13:23:57	29/01/95

The measuring electrode should be connected to the coaxial INPUT connector. Its banana plugs should be inserted in the °C terminals.

## *Inputs*

# MEASUREMENTS

### **Principle:**

## *Theory*

The oxygenmeter and its electrode function according to the Clark principle with platinum as cathode and silver as anode in an electrolyte cell to which a polarisation voltage of 800 mV is applied. Oxygen gas present in the electrolyte is reduced to OH ions at the platinum cathode. The resulting current is diffusion limited and therefore proportional to the oxygen concentration in the sample solution. This current is amplified, corrected, and displayed in mg/l, ppm or % dissolved oxygen.

### **Electrode:**

The anode and cathode of the electrode are immersed in a special buffer solution inside the electrode, to give high stability readings and short polarisation times. The electrolyte compartment is separated by a PTFE film from the sample solution. Only gases, but no ions, can diffuse through this film from the sample solution into the internal electrolyte.

### **Temperature compensation:**

Speed of electrode reactions depends on temperature. Therefore, for a given oxygen concentration, the electrode current becomes greater at higher temperature, and conversely. This temperature effect is compensated with the NTC resistor incorporated in the electrode. For aqueous solutions the oxygen meter therefore always gives correct readings in ppm, independent of the actual temperature.

## Interferences:

All substances which can diffuse through the membrane and for which 800 mV potential suffices for polarographic reduction, will be reduced in the electrode. This will give a corresponding current contribution, if they are present. Interference can be caused by ions entering the electrode through porous or mechanically damaged membranes and by diffusion of other reactive gases apart from oxygen, e.g. CO<sub>2</sub>, Cl<sub>2</sub>, SO<sub>2</sub>, and H<sub>2</sub>S. These substances react in undesired manner with the electrode.

Acidic or basic gases change the pH value of the electrolyte solution and thus disturb the reading, particularly when measuring small oxygen concentrations. High salt concentrations in the sample solution can falsify readings too.

Always store the electrode in distilled water after use! The membrane may cease to function correctly after prolonged use or mechanical damage. This will falsify the readings and the electrolyte may drain out. To replace the membrane (see page 49), unscrew the measuring head and dismantle it further to remove the old membrane. Take a new membrane and mount it carefully, fill the measuring head with electrolyte solution, and then slowly screw it onto the electrode. The electrode is now ready for use again.

The electrode body consisting of the electrode shaft and the plug head has almost unlimited life and requires no maintenance. After prolonged use of the electrode in solutions containing high concentrations of oxygen, the silver anode may become deactivated by undesired deposits. An indication is that the electrode no longer responds correctly to calibration. In this case, screw off the measuring head and cautiously remove the precipitated deposits from the silver electrode surface with the aid of filter paper. The electrode is then ready again for operation after refilling the measurement head with new electrolyte solution.

## Start this method only with a well polarised electrode!

1. Before making measurements, check that the membrane is undamaged and that there are no gas bubbles present in the electrolyte space.
2. Connect the electrode and switch the instrument on. After a few minutes, the electrode is polarised and ready for making measurements.
3. Select the ppm (mg/l) or % range by pressing **MODE**. The display will immediately show the measured oxygen according to the previous calibration. Should you want to restandardise, press **CAL**.
4. The display shows the salinity correction, e.g. [SAL=2.1%]. Leave this value to zero unless you are going to measure in heavily salted solutions such as e.g. sea-water (3.5%). Pre-set the proper value with ▲ or ▼ and press **CAL** or **MODE** to cancel.
5. The display shows the pressure correction, e.g. [P=1023hP]. Pre-set the actual air pressure (hP) with ▲ or ▼ and press **CAL** or **MODE** to cancel.
6. Press ▲ or ▼ until the display shows [1.Point]. Press **CAL** to start the calibration procedure.

## ***Maintenance***

## ***Measurements with 1-point calibration***

7. The electrode exposed in this manner to the air reaches an equilibrium corresponding to the partial pressure of oxygen and thus to saturation in water at the given temperature. Alternatively, rinse the electrode with distilled water and then dip it into freshly air saturated distilled water for calibration. The advection rate must be at least 10 cm/s (**stir!**). Air saturated water can be prepared by shaking or vigorously stirring distilled water for 10 minutes with free access of air. The instrument shows the measured saturation, e.g. [SA<sub>t</sub>=8.3] and will calibrate automatically when readings are stable ([=] stops blinking).
8. After rinsing the electrodes with distilled water, immerse them in the samples and read the display. The advection rate must be at least 10 cm/s (**stir!**).

1. Before making measurements, check that the membrane is undamaged and that there are no gas bubbles present in the electrolyte space.
2. Connect the electrode and switch the instrument on. After a few minutes, the electrode is polarised and ready for making measurements.
3. Select the ppm (mg/l) or % range by pressing **MODE**. The display will immediately show the measured oxygen according to the previous calibration. Should you want to restandardise, press **CAL**.
4. The display shows the salinity correction, e.g. [SAL=2.1%]. Leave this value to zero unless you are going to measure in heavily salted solutions such as e.g. sea-water (3.5%). Pre-set the proper value with ▲ or ▼ and press **CAL** or **MODE** to cancel.
5. The display shows the pressure correction, e.g. [P=1023hP]. Pre-set the actual air pressure (hP) with ▲ or ▼ and press **CAL** or **MODE** to cancel.
6. Press ▲ or ▼ until the display shows [2.Point]. Press **CAL** to start the calibration procedure.
7. Dip the electrode into a saturated sodium sulphite solution (without stirring). The instrument shows the measured zero-current (nA) of the electrode, e.g. [O<sub>2.o</sub>=1.3] and will calibrate automatically when readings are stable ([=] stops blinking).
8. Rinse the electrode with distilled water and cautiously dry it without rubbing (use cotton wool or soft filter paper). The electrode exposed in this manner to the air reaches an equilibrium corresponding to the partial pressure of oxygen and thus to saturation in water at the given temperature. Alternatively, rinse the electrode with distilled water and then dip it into freshly air saturated distilled water for calibration. The advection rate must be at least 10 cm/s (**stir!**). Air saturated water can be prepared by shaking or vigorously stirring distilled water for 10 minutes with free access of air. The instrument shows the measured saturation, e.g. [SA<sub>t</sub>=8.3] and will calibrate automatically when readings are stable ([=] stops blinking).
9. After rinsing the electrodes with distilled water, immerse them in the samples and read the display. The advection rate must be at least 10 cm/s (**stir!**).

### ***Measurements with 2-point calibration***

### Relative O<sub>2</sub> measurement:

1. While in the O<sub>2</sub> range, press  $\blacktriangle$ . At that moment the instrument stores the actual value and automatically subtracts it from all next measurements. You can repeat this as many times as required. A blinking [r] in the display appears.
2. To cancel relative readings, press **MODE** and you will be in the normal measuring mode again.

- \* You can also check the saturation-current (nA) and the saturation-point (ppm) by pressing  $\blacktriangledown$  sequentially: e.g. [Cur=67.9] & [Sat=8.73]. See also GLP-protocol.
- \* The instrument will refuse automatic standardisation when the electrode is unstable. A worn electrode may be the cause. Select a lower resolution for a more rapid standardisation or replace the membrane and electrolyte of the electrode!
- \* After replacement or refilling of the measuring head the first polarisation of the electrode could last  $\pm 15$  minutes!
- \* Press **RES** to change the resolution from 0.1 to 0.01 ppm or from 1 to 0.1 %.

### *Relative measurement*

### Measurement:

1. Read temperature by pressing  $^{\circ}\text{C}$ .

### *Temperature measurement*

### Calibration:

1. Press **CAL** to calibrate the  $^{\circ}\text{C}$  scale with any reference. Immerse the electrode in a solution of known temperature.
2. The display shows e.g. [ $^{\circ}\text{C}=22.3$ ]. Calibrate to the proper value with  $\blacktriangle$  or  $\blacktriangledown$  and press **CAL** again or press **RES** to reset the absolute readings.

## DATA COMMUNICATION

A recorder can be connected to the (+) and (-) terminals. The 0 to 2 V output range corresponds with 0 to 60 ppm or 600 %. Use only laboratory recorders with a high input impedance!

### *Recorder output*

1. Select the O<sub>2</sub> range (ppm or %) by pressing **MODE**.
2. Check the saturation-current (nA) and the saturation-point (ppm) by pressing  $\blacktriangledown$  sequentially: e.g. [Cur=67.9] & [Sat=8.73].
3. Press **PRINT** to print a complete calibration report, e.g.:

### *GLP-protocol*

CONSORT Oxygenmeter Z621

-----  
SETTINGS

-----  
Identification No : 004  
Software Version : 3.0  
User Code : off  
Relative O<sub>2</sub> : off

NTC Cal.Temp. (C) : 30.0

Salinity (%) : 0.0  
Pressure.(hP) : 1005

Alarm O<sub>2</sub> : on  
Alarm High (ppm) : 10.7  
Alarm Low (ppm) : 5.0

Batt.Capacity (%) : 87

CALIBRATION

-----  
Cal. points : 2

Salinity (%) : 0.0  
Pressure (hP) : 1035  
Temperature (C) : 19.4  
Saturation (ppm) : 7.48

Zero-point (nA) : 1.2  
Response Time (s) : 96

STATISTICS

-----  
Salinity (%) : -0.3  
Pressure (hP) : +25  
Temperature (C) : -1.4

Zero-point (nA) : +0.2  
Response Time (s) : -16

A standard RS232 output terminal (DP9) is provided for interfacing the instrument with a printer or computer. Data is sent in the ASCII code at a BAUD rate of 2400 b/s (8 bit, no parity, 1 start and 2 stopbits, automatic line feed ON, data transfer protocol Xon/Xoff).

***Digital output***

### Serial port pinout specifications:

pin 1: RLSD, received line signal detector  
pin 2: TxD, transmit data  
pin 3: RxD, receive data  
pin 4: not connected  
pin 5: Gnd, signal ground  
pin 6: DSR, data set ready  
pin 7: not connected  
pin 8: CTS, clear to send  
pin 9: not connected

### Interfacing with a printer:

Press **PRINT** each time you want to print the measured values. Run the 'printing data' procedure (page 6) for automatic printing at timed intervals, while you are measuring. The **PRINT** keypath does not function in this case. It is also disabled during an automatic data-logging.

### **Printer connection**

#	VALUE	UNIT	C
0001	7.34	ppm	23.7
0002	7.21	ppm	23.7
0003	7.08	ppm	23.8
0004	6.71	ppm	23.8
0005	6.49	ppm	23.9

### Using WINDOWS™ to collect data:

1. Connect the computer with the RS232 terminal, using the optional AK2310 cable (eventually with adapter AB3190 for connection with a 25 pole RS232 terminal).
2. Start **WINDOWS™** and select *Terminal* in the *Accessories Group*.
3. Choose *Receive Text File* in the *Transfer Menu* and create a file name (TXT format) for the data to receive.
4. Select the *Settings Menu* and make sure the following parameters are programmed:

Baud Rate : 2400  
Data bits : 8  
Stop bits : 2  
Parity : None  
Flow control: Xon/Xoff

5. Start the oxygenmeter in the desired measuring mode for manual or automatic data collection (see 'Data acquisition' page 4 for more details).
6. When finished, open the created file with a standard spreadsheet (e.g. **EXCEL™**) or wordprocessor (e.g. **WORD™**) for further processing.

### **Computer connection**

## Interfacing with a computer:

Send a single character to the instrument and it will execute the corresponding command as follows:

- "1" = press **MODE**.
- "2" = press °C.
- "3" = press ▲.
- "4" = press ▼.
- "5" = press **CAL**.
- "6" = press **HOLD**.
- "7" = press **RES**.
- "8" = press **PRINT**.
- "?" = send display to computer.
- "+" = keyboard on (manual operation remains possible).
- "-" = keyboard off (manual operation disabled).

Follow exactly the same measuring or calibration procedures, as described in this manual, to include any desired command in your computer data-acquisition program.

This example continuously prints the measured values on the computer screen:

10 ON ERROR GOTO 70	'makes sure no error occurs
20 OPEN "COM1:2400,N,8,2" AS #1	'prepares computer input
30 PRINT #1,"8"	' <b>PRINT</b> button pressed
40 INPUT #1,V\$	'measured values transmitted
50 PRINT V\$	'received values printed
60 GOTO 30	'repeat procedure
70 RUN	

# TROUBLE SHOOTING

*This procedure permits to calibrate automatically all amplifiers of the instrument.*

## Calibrate

1. Leave all inputs open.
2. Switch the instrument ON while holding **CAL** pressed.
3. The display shows [SELFCAL] for a while and returns to the measurements.

*This procedure resets the instrument to the original factory settings.*

## Reset

1. Switch the instrument ON while holding **MODE** pressed.
2. The display shows [rESEt] for a while and returns to the measurements.

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