C32xx

MANUAL HANDLEIDING MODE D'EMPLOI ANLEITUNG

Rev. 04/2020

C3210
C3230

C3210

CH.1: pH, mV, Conductivity + Ref. + °C
CH.2: pH, mV, Conductivity + Ref. + °C

C3230

CH.1: pH, mV, Ion, O2, Conductivity + Ref. + °C
CH.2: pH, mV, Ion, O2, Conductivity + Ref. + °C

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General Information

Introduction

This instrument is manufactured with the latest technology and needs no particular maintenance. **Consort** certifies that this instrument was thoroughly inspected and tested at the factory prior to shipment and found to meet all requirements defined by contract under which it is furnished. However, dimensions and other physical characteristics may differ.

The normal operating temperature should be between 0° and 40°C. Never use the instrument in a room with high humidity (>95 %) or at very low temperatures (condensation water!).

Manufacturer

Consort BV
Hertenstraat 56/9 Tel (+32)(14)41 12 79
B2300 Turnhout Fax (+32)(14)42 91 79
Belgium E-mail: info@consort.be

Warranty

This instrument (excluding all accessories) is warranted against defective material and workmanship for period of thirty-six (36) months from the date of shipment ex factory. **Consort** will repair all defective equipment returned to it during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures in this instruction manual and when not having been subject to accident, alteration, misuse or abuse. A return authorisation must be obtained from **Consort** before returning any product for warranty repair on a freight prepaid basis!

Consort is not liable for consequential damages arising out of the use or handling of its products.

Servicing

In the event of this instrument being returned for servicing, the owner is requested **NOT** to send the following items unless they are suspect:

Manual
Cables
Accessories

If serious malfunctioning occurs, stop using the unit immediately and consult your local **Consort** dealer.

AC adaptor

Connect the mains adapter to the DC socket. Do not hold the adapter by wet hand.

Keyboard

MODE	=	Selects the settings or escapes from error traps, calibration procedures, etc..by returning to the original mode.
↑↓ selecting a function	=	Buttons for browsing between the selected modes, for entering a value or for value
OK	=	Proceeds a function.
CAL	=	Starts a calibration.
STORE	=	Stores the displayed value or send it to a computer or printer.
HOLD	=	Holds display when measuring.

Installation of the instrument

Inputs

The measuring electrodes should be connected to the coaxial connectors CH1 or CH2. Automatic temperature compensation and measurement are possible by connecting a Pt1000 temperature compensator to the banana connectors °C1 or °C2. Without compensator, the manual temperature compensation is automatically switched on.

Outputs

A standard USB port is provided for interfacing the instrument with a computer or network. A standard RS232 port is provided for interfacing the instrument with a printer. Data is sent in the ASCII code at a BAUD rate of 1200...115200 bps (8 bit, no parity, 1 stopbit).

Language

1. Press **MODE**.
2. Select [SYSTEM] and press **OK**.
3. Select [LANGUAGE] and press **OK**.
4. Select the desired language and press **OK**.

Contrast

1. Press **MODE**.
2. Select [SYSTEM] and press **OK**.
3. Select [CONTRAST] and press **OK**.
4. Adjust the contrast and press **OK**.

Date and time

1. Press **MODE**.
2. Select [SYSTEM] and press **OK**.
3. Select [DATE/TIME] and press **OK**.
4. Select [MODIFY] and press **OK**.
5. Adjust date and time and press **OK**.

Password

1. Press **MODE**.
2. Select [SYSTEM] and press **OK**.
3. Select [PASSWORD] and press **OK**.
4. A private code can be programmed to avoid undesired access to the instrument. Select [YES] and press **OK** to enter your secret sequence of 5 keys.

Stand by

1. Press **MODE**.
2. Select [STAND BY] and press **OK**.
3. Select [YES] (= continue automatically after a power failure) or [NO] (= instrument can be switched off).

Channels

1. Press **MODE**.
2. Select [CHANNELS] and press **OK**.
3. Select the desired channel and press **OK**.
4. Select the desired range and press **OK**.
5. Select the desired input for temperature compensation and press **OK**.

User standard tables

- Press **MODE**.
- Select [USER TABLES] and press **OK**.
- Select [pH buffers] or [S/cm standards] and press **OK**.
- Follow the instructions on the screen to enter the different values. The value for 25°C is obligatory!

pH/mV/temperature measurement

pH measurement

- Select the desired mode (pH) by pressing $\uparrow\downarrow$. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
- The instrument permits to choose between the buffers in memory (1.68, 2.00, 4.00, 4.01, 6.87, 7.00, 9.18, 9.21, 10.01, 12.00, 12.45 + up to 5 user tables). Select the proper values and press **OK**. The unused buffers should be switched off.
- Rinse the electrodes with distilled water and immerse them in the first buffer solution. Select [CALIBRATE], press **OK** and follow the instructions on the screen until the calibration is finished.
- After rinsing the electrodes with distilled water, immerse them in the samples and read the display.
- Rinse the electrodes always with distilled water after use and store them in a 3...4 M KCl solution.
- *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
- *Stirring the solution during the measurements promotes the homogeneity and is obligatory!*
- *The instrument will refuse automatic calibration when the electrode is unstable. Insufficient stirring or a worn electrode may be the cause.*

Using pH electrodes with an abnormal zero point

1. Press **CAL**.
2. Select [Eo = x mV] and press **OK**.
3. Enter the zero point of the electrode and press **OK**.
4. *Hold MODE pressed and press \downarrow to change the resolution from 0.1 to 0.001 pH.*

mV measurement

1. Select the desired mode (mV) by pressing $\uparrow\downarrow$. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
2. Immerse the electrodes in a standard solution of known potential. Select [CALIBRATE] and press **OK**. Adjust to the proper value and press **OK**.
6. *Select [RESET] and press **OK** to reset the calibration.*
7. *Hold MODE pressed and press \hat{e} to change the resolution from 1 to 0.1 mV.*

Temperature measurement

1. Select the desired mode ($^{\circ}\text{C}$) by pressing $\hat{e}\hat{e}$. Should you want to recalibrate, press **CAL**.
2. Immerse the Pt1000 in a solution of known temperature. Select [CALIBRATE] and press **OK**. Adjust to the proper value and press **OK**.
- *Select [RESET] and press **OK** to reset the calibration.*
- *Without Pt1000, press **CAL**, adjust the manual temperature compensation and proceed by pressing **OK**.*

Conductivity/TDS/Salinity/Resistivity measurement

Conductivity measurement

1. Select the desired mode (S/cm) by pressing $\uparrow\downarrow$. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
 2. The instrument permits to choose between the standards in memory (1413 $\mu\text{S/cm}$, 12.88 mS/cm, 111.8 mS/cm + up to 3 user tables). Select the proper values and press **OK**. The unused standards should be switched off.
 3. After rinsing the electrode several times with the first standard solution, immerse it in that standard. The solution temperature is not so critical but should lie between 0°C and 40°C. When no Pt1000 is used, do not forget to compensate manually first! Select [CALIBRATE], press **OK** and follow the instructions on the screen until the calibration is finished.
 4. Rinse the electrode several times with the sample, immerse it in that solution and read the display.
 5. Rinse the electrode always after use and store it in distilled water (add some detergent to keep the spongy platinum surface in perfect condition).
1. *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
 2. *Stirring the solution during the measurements promotes the homogeneity and is obligatory!*
 3. *The instrument will refuse automatic calibration when the electrode is unstable. Insufficient stirring or a worn electrode may be the cause.*
 4. *When starting a titration, hold **MODE** pressed and press \hat{e} to lock the actual range and avoid cross-over errors due to differences in the measuring frequency. Hold **MODE** pressed and press \downarrow to return again to the normal measuring mode.*

Temperature compensation

1. Press **CAL**.
2. Select [REF.] and press **OK**.
3. Select the temperature to which all future measurements will be referred to (25°C or 20°C) (C305x: 20°C...40°C in 1°C steps) and press **OK**. Select the value 0 to disable the temperature compensation and read absolute conductivity measurement values.

Capacitive compensation

1. Capacitive compensation increases accuracy in the very low conductivity ranges (<10 $\mu\text{S/cm}$). Select whether or not this compensation should be applied. Verify if the attached electrode is completely dry and press **CAL**.
2. Select [CAP.COMP] and press **OK**.
3. Select [YES] or [NO] and press **OK**. Follow the instructions on the screen.

TDS measurement

1. Select the desired mode (mg/l)(TDS) by pressing $\uparrow\downarrow$. The display will immediately show the measured value.

Salinity measurement

1. Select the desired mode (SAL) by pressing $\uparrow\downarrow$. The display will immediately show the measured value.

Resistivity measurement

1. Select the desired mode (W.cm) by pressing $\uparrow\downarrow$. The display will immediately show the measured value.
- To browse between conductivity, resistivity, salinity or TDS, hold **MODE** pressed and press $\square\square$ sequentially.

Dissolved oxygen/Air pressure measurement

Dissolved oxygen measurement

1. Select the desired mode (ppm) or (%O₂) by pressing **↑↓**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
2. The display shows the salinity correction. Select the proper value and press **OK**. Leave salinity correction to zero unless you are going to measure in heavily salted solutions such as e.g. sea-water (35 g/l). Select [CALIBRATE], press **OK** and follow the instructions on the screen until the calibration is finished.
3. The electrode exposed to the air reaches an equilibrium corresponding to the partial pressure of oxygen and thus to saturation in water at the given temperature. The instrument shows the measured saturation, current, temperature, and will calibrate automatically when readings are stable.
4. After rinsing the electrodes with distilled water, immerse them in the samples and read the display. Stirring the solution during the measurements promotes the homogeneity and is **obligatory!** The advection rate must be at least 10 cm/s.
5. Rinse the electrode always after use and store it in distilled water.
 - *A blinking decimal point warns you for unstable measurements. Wait to read the display!*
 - *Stirring the solution during the measurements promotes the homogeneity and is therefore always recommended.*
 - *Hold **MODE** pressed and press **↓** to change the resolution from 0.1 to 0.01 ppm or from 1 to 0.1 %.*

Air pressure measurement

1. Select the desired mode (hPa) by pressing **↑↓**. Should you want to recalibrate, press **CAL**.
2. Select [CALIBRATE], press **CAL**, adjust to the real air pressure and press **OK**.
 - *Select [RESET] and press **OK** to reset the calibration.*

Principle

The oxygen meter and its electrode function according to the Clark principle with silver as cathode and lead as anode in an electrolyte cell. Oxygen gas present in the electrolyte is reduced to OH ions at the cathode. The resulting current is diffusion limited and therefore proportional to the oxygen concentration in the sample solution. This current is amplified, corrected, and displayed in mg/l, ppm or % dissolved oxygen.

Interferences

All substances which can diffuse through the membrane and for which 800 mV potential suffices for polarographic reduction, will be reduced in the electrode. This will give a corresponding current contribution, if they are present. Interference can be caused by ions entering the electrode through porous or mechanically damaged membranes and by diffusion of other reactive gases apart from oxygen, e.g. CO₂, Cl₂, SO₂, and H₂S. These substances react in undesired manner with the electrode. Acidic or basic gases change the pH value of the electrolyte solution and thus disturb the reading, particularly when measuring small oxygen concentrations. High salt concentrations in the sample solution can falsify readings too.

Ion selective measurement (C3230 only)

Calibration and measurement

1. Select the desired mode by pressing **↑↓**. The display will immediately show the measured value according to the previous calibration. Should you want to recalibrate, press **CAL**.
2. The instrument permits to choose between the possible standards (1, 10, 100 ng/l, 1, 10, 100 µg/l, 1, 10, 100 mg/l, 1, 10, 100 g/l). Select the proper values and press **OK**. The unused standards should be switched off.
3. Select [CALIBRATE] and press **OK**.
4. After rinsing the electrodes with distilled water, immerse them in the first standard solution. Select the first calibration value and press **OK** when readings are stable. Go on in the same way with the next standard solutions.
5. Decide whether a blank correction should be carried out or not. Follow the instructions on the screen. After rinsing the electrodes with distilled water, immerse them in a blank solution. When readings are stable press **OK**.
6. Rinse the electrodes with distilled water, immerse them in the samples, and read the concentration on the display.
5. *Maintenance: we refer to the manual supplied with the ion specific electrodes.*

Good Laboratory Practice

1. Select the desired range by pressing **↑↓** and then press **CAL**.
2. Select [GLP] and press **OK**.
3. Select [SHOW REPORT] and press **OK**. Browse with **↑↓** to show a complete calibration report.
4. Select [SEND REPORT] and press **OK** to send the report to a computer.

Calibration reminder

1. Select the desired range by pressing **↑↓** and then press **CAL**.
2. Select [GLP] and press **OK**.
3. Select [INTERVAL] and press **OK**.
4. Select the desired time interval between each automatic warning for a new calibration of the electrodes and press **OK**.

Data Storage in the internal memory

1. Press **MODE**.
2. Select [DATA] and press **OK**.

Start the data-logging

1. Select [ACTIVATE] and press **OK**.
2. Decide whether data-logging should be carried out or not and press **OK**.
3. Select [CONTINUOUS] and press **OK**.
4. Decide whether the oldest data should be overwritten when the memory limit is reached or not and press **OK**.
5. Select [INTERVAL] and press **OK**.
6. Select the desired time interval between the data-logging and press **OK**.
7. Return to the measurements by pressing **MODE**.
8. The data-logging starts according to the previous settings while the display shows the logging-number e.g. [LOG.00027].

Process the stored values

1. Select [PROCESS] and press **OK**.
2. Select [TABLE] or [SEND] to display or send the stored data and press **OK** to continue. Follow the instructions on the screen.

Erase the stored values

1. Select [PROCESS] and press **OK**.
2. Select [ERASE] and press **OK**. Follow the instructions on the screen.

External storage in a computer using USB or RS232

1. Press **MODE**.
2. Select [SERIAL PORT] (only on Ethernet versions) and press **OK**.
3. Select [INTERVAL] and press **OK**.
4. Select the desired interval between the transmitted data and press **OK**. Pre-set to zero if no automatic transmitting is required.
5. Select [BAUD] and press **OK**.
6. Select the desired transmission rate and press **OK**.
- *You can download free software and a suitable USB driver or an Ethernet configuration document from www.consort.be.*

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DECLARATION OF CONFORMITY

We declare under our sole responsibility that the product

**Multi-channel analysers
content of the type numbers**

C3210, C3220

to which this declaration relates is in conformity
with the following standards

LOW VOLTAGE DIRECTIVE 2006/95/EC

EN61010-1

EMC DIRECTIVE 2004/108/EC

EN61326-1

Consort BV

Hertenstraat 56/9, B-2300 Turnhout, Belgium

Tel: (+32) (0)14 41 12 79

Fax: (+32) (0)14 42 91 79

Sales: sales@consort.be

The electro-magnetic susceptibility has been chosen at a level that gains proper operation in residential areas, on business and light industrial premises and on small-scale enterprises, inside as well as outside of buildings. All places of operation are characterised by their connection to the public low voltage power supply system.

WARRANTY CERTIFICATE

This instrument (excluding all accessories) is warranted against defective material and workmanship for a period of thirty-six (36) months from the date of shipment ex factory.

Accessories and breakable items such as electrodes are not warranted unless proven to be defective before shipment.

The original purchase order numbers, Consort invoice numbers and serial numbers of the products must be provided.

CONSORT will repair all defective equipment returned to it during the warranty period without charge (CIF Turnhout prepaid by sender), provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures described in the instruction manual and when not having been subject to accident, alteration, misuse or abuse.

If the products have been used with or have come into contact with fluids, an MSDS (material safety data sheet) must be supplied prior to issuing a return authorisation.

A return authorisation must be obtained from **CONSORT** before returning any product for warranty repair on a freight prepaid basis!

Consort BV

Hertenstraat 56/9
B-2300 Turnhout • Belgium

Tel : (+32) (0)14 41 12 79
Fax : (+32) (0)14 42 91 79

Sales : sales@consort.be
Support : support@consort.be
Information : info@consort.be